

Energy Transfer - Wind Turbine

What's included


- Fan
- Spacer
- · Thumb nut

Equipment setup

The Wind Turbine accessory is designed to be attached to the Energy Transfer Generator to create a wind energy turbine. There are two versions of the Energy Transfer Generator: ET-8771A and ET-8771B. Follow the directions below for your version of the Energy Transfer generator.


ET-8771B instructions

- 1. Remove the knob and three-step pulley from the ET-Generator.
- 2. Slide the spacer that comes with the Wind Turbine onto the generator shaft.
- Push the Wind Turbine onto the generator shaft so that the turbine blades curve away from the generator. Note that the Wind Turbine fits tightly on the generator shaft.
- 4. Screw the ET-Generator knob onto the shaft. Do not use the thumb nut that comes with the Wind Turbine.

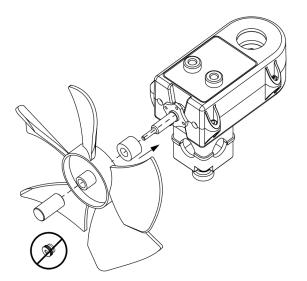

ET-8771A instructions

- Remove the thumbnut and three-step pulley from the ET-Generator.
- 2. Slide the spacer that comes with the Wind Turbine onto the generator shaft.
- Push the Wind Turbine onto the generator shaft so that the turbine blades curve away from the generator. Note that the wind turbine fits tightly on the generator shaft.
- Screw the thumbnut that comes with the Wind Turbine onto the shaft.

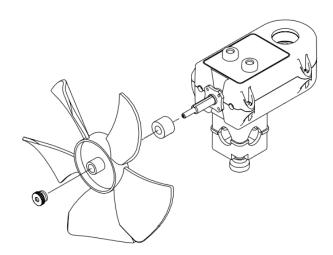

Technical Support

Need more help? Our knowledgeable and friendly Technical Support staff is ready to answer your questions or walk you through any issues.

Chat pasco.com

Phone 1-800-772-8700 x1004 (USA)

+1 916 462 8384 (outside USA)

Warranty

For a description of the product warranty, see the Warranty and Returns page at pasco.com/legal.

Trademarks

PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information, visit pasco.com/legal.