# **PASCO**Collision Cart

## MF-9454

The Collision Cart is a 500-gram cart with an aluminum body and ABS end caps. It is the companion to the Plunger Cart (ME-9430), which is blue and has a spring plunger. The different colors make it easier to keep track of each cart during collision studies. The cart includes magnets, and hook and loop pads for exploring elastic and inelastic collisions. The cart is compatible with all PASCO Dynamics Tracks and accessories.

## Part descriptions



#### 1. Wheels

The wheels are mounted to low-friction ball-bearings. The wheels retract into the body so that they are protected from drops or being stepped on. The wheels are designed to fit in the grooves on PASCO tracks.

## 2. ABS End Caps

## 3. Magnets

Use for elastic collision experiments.

## 4. Upper Tie Point

Use for attaching a string or springs (ME-8999) to the cart.

#### Slot

Use for inserting a Cart Picket Fence (ME-9804).

## 6. Mounting Holes

Use the threaded holes to attach various accessories.

## 7. Accessory Tray

Use to hold Cart Masses (ME-6757A).

## 8. Hook and Loop Pads

Use for inelastic collision experiments.

## 9. Lower Tie Point

Use for attaching a string to pull the cart.

# Replacing the axles

The bearings inside the axles may become damaged from dirt and other contaminants over time. Replacement axles are available as item ME-6957.

## Additional resources

## **Product information**

Visit the product web page at pasco.com/product/ME-9454 for additional information including:

- Specifications
- Buying Guide
- Experiments
- Documents



### **Experiments**

Print-ready experiment worksheets are available to download from the PASCO website. Go to pasco.com/freelabs and enter ME-9454 in the Part No. field.

## **Technical Support**

Need more help? Our knowledgeable and friendly Technical Support staff is ready to provide assistance with this or any other PASCO product.

Phone (USA) 1-800-772-8700 (Option 4)

Phone (International) +1 916 462 8384

Online pasco.com/support

# Warranty, Copyright, and Trademarks

## Limited Warranty

For a description of the product warranty, see the Warranty and Returns page at www.pasco.com/legal.

#### Copyright

This document is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

### Trademarks

PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.