

OS-7054; Meter Stick Optics Battery Replacement Instructions

Battery part number PS-3295; Li-Poly 1000 mAh

Recommended Equipment: (PS-3299; Battery Replacement Tool Set)



Torx T8 Driver




Torx T5 Driver



PS-3295




CAUTION: Failure to use proper ESD protection could result in damage to the electronics.

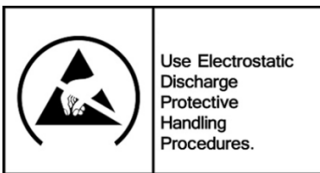


Step 1: Position the Device on a surface as shown. Note the location of the (4) Torx T8 Screws.

Torx T8 Screw, 4 Places

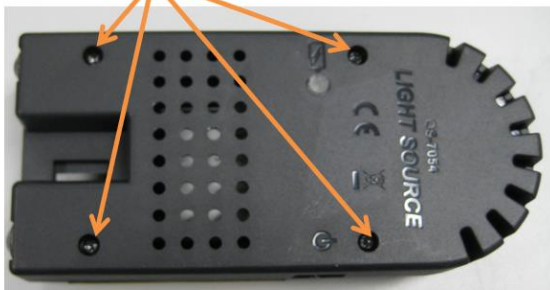


Step 2: Unscrew and remove the (4) Screws with the Torx T8 Screwdriver.

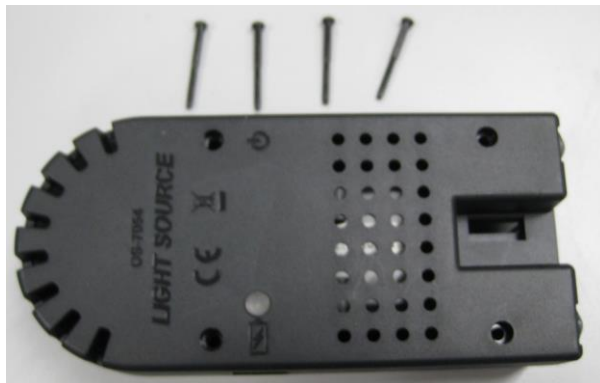


PASCO Form Number	Description	Date		Revision
013-17796	OS-7054 Battery Replacement Instructions	4-30-24	Page 1 of 3	A

Step 3 Remove the back and set aside. Note the location of the (2) Torx T5 Screws holding the PC Board. Remove the (2) screws with a Torx T5 Screwdriver and lift the PC Board as shown.

Torx T5 Screw, 2 Places

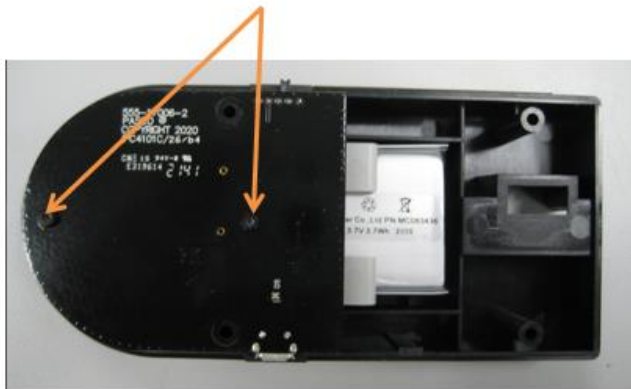

Step 4: Unplug the battery from its connector. Lift off the plastic hood and set aside.

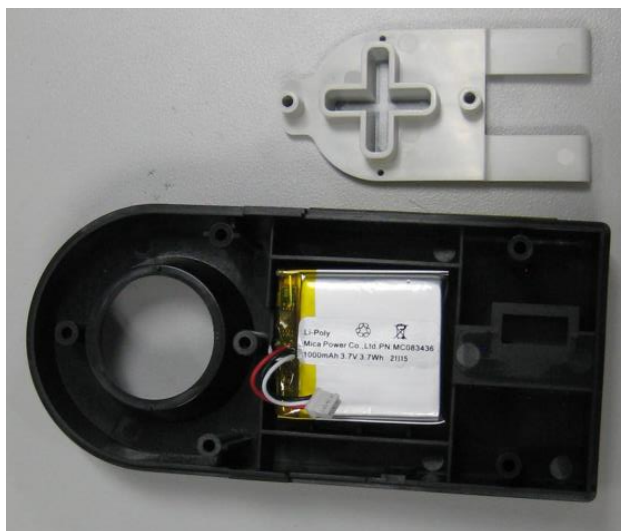

Step 5: Remove the battery and add the replacement battery. Plug the battery into its mating connector and reverse the steps noted above to complete the reassembly.

PASCO Form Number	Description	Date		Revision
013-17796	OS-7054 Battery Replacement Instructions	4-30-24	Page 2 of 3	A

Battery Replacement and Disposal Instructions (Reference):



Batteries contain chemicals that, if released, may affect the environment and human health. Batteries should be collected separately for recycling, and recycled at a local hazardous material disposal location adhering to your country and local government regulations. To find a battery recycling drop-off location, please contact local waste disposal service or the product representative.



The battery or batteries used in PASCO products are marked with the European Union symbol for waste batteries (above) to indicate the need for separate collection and recycling. For small batteries, the symbol is printed on the packaging.

If you need further assistance, please contact Technical Support at support@pasco.com or through www.pasco.com on chat. To access chat through your browser, please enable all cookies on our website.

PASCO Form Number	Description	Date		Revision
013-17796	OS-7054 Battery Replacement Instructions	4-30-24	Page 3 of 3	A