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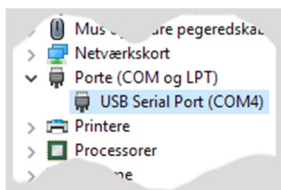
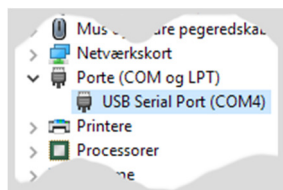
Frederiksen's Geiger-Müller counter 513610 (as well as older version 513600) can be used with the free software *Datalyse*. The communication goes through a 512565 USB communication adapter for 513610.

512565 is a USB to serial adapter. Internally in the PC, it will appear as a serial port a.k.a. a COM port. When installed correctly, the adapter will appear in Windows *Device Manager* under *Ports* when it is plugged in.

On the image you can see that the port in this case is called COM4.

If instead the adapter appears under *Unknown devices*, e.g. with the name "TTL-232R-3V3-AJ", a driver must be installed for the device.

In Windows 10, you can simply double-click at the line with the device and select *Update Driver*. In the following dialog you choose to install the driver automatically (the PC must have Internet access).

Older versions of Windows may require a different procedure – we are not covering that here.

Download Datalyse

The program is free and can be downloaded here:

www.datalyse.dk

It exists in an English version. The program doesn't need any installation. You start it by double-clicking the program file.

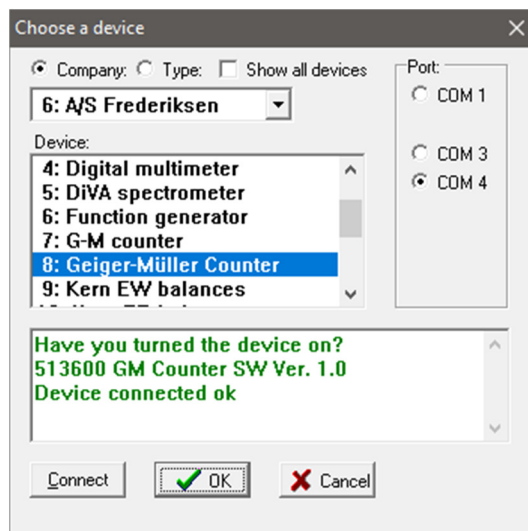
Select device and port in Datalyse

Select in the menu: *Device / Choose device*.

In the drop-down menu for companies, select "A/S Frederiksen"

On the *Device* list, choose "Geiger-Müller Counter"

There may be more ports to choose among – pick the one you found in *Device Manager*.

Now, the program is ready to receive data from the Geiger counter.

Support

Frederiksen Scientific recommends the use of *Datalyse* with e.g. our Geiger counters. It is a very useful program and easy to operate. But:

Frederiksen Scientific did not develop the *Datalyse* software and it is not sold by us. We can therefore not offer end-user support for using it.

Documentation can be found on the homepage for *Datalyse*.

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I dropdown-menuen for firmaer vælges "A/S Frederiksen"