

## Quick Start

### What You Need

ScienceWorkshop 750  
USB Interface



USB A-B cable



AC adapter



DataStudio CD  
(version 1.7 or later)




### Install Software

- 1 Insert the DataStudio CD into your computer's CD-ROM drive.
- 2 The DataStudio Installer will automatically begin the installation process.
- 3 Follow the on-screen prompts.



### Get Connected

- 1 Connect the power supply to the 750 interface and turn on the interface.
- 2 Insert the "B" end of the USB cable into the USB port on the back of the interface (see below).
- 3 Insert the "A" end of the USB cable into a USB port on the computer (see below).
- 4 Open DataStudio to begin an experiment.

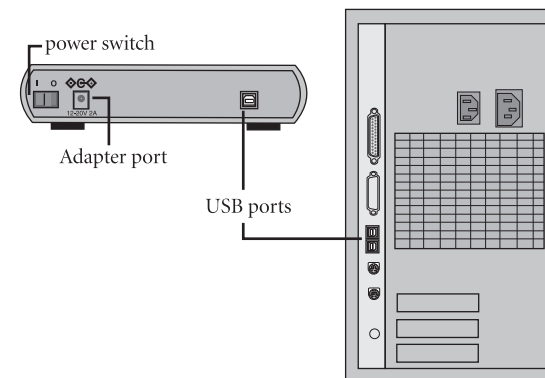

For free technical assistance, please call our technical support department

fax **1.916.786.3292**

012-08164A

email **techsupp@pasco.com**

Internet **www.pasco.com**

Telephone 7:00am - 4:30pm PST

In the U.S. **800.772.8700**

International **1.916.786.3800**