

Quick Start

What You Need

850 Universal Interface



USB
cable



power supply
and
power cord



PASCO Capstone software




Install Software

- 1 Insert the PASCO Capstone software storage device into your computer.
- 2 The Capstone Installer will automatically begin the installation process.
- 3 Follow the on-screen prompts.

Get Connected

- 1 Connect the power supply to the power port on the back of the 850 interface and turn on the interface.
- 2 Insert the small end of the USB cable into the mini USB port on the back of the interface (see below).
- 3 Insert the larger end of the USB cable into a USB port on the computer (see below).
- 4 Open Capstone to begin an experiment.

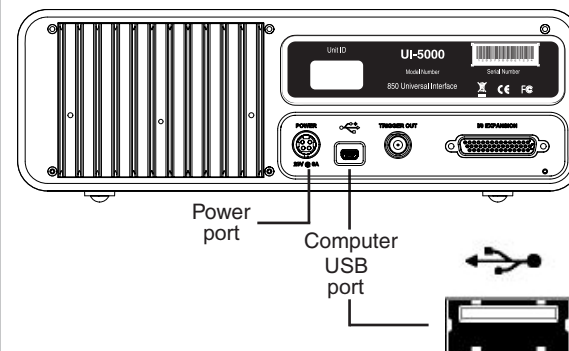

For free technical assistance, please call our technical support department

fax **1.916.786.7565**

012-13507A

email **techsupp@pasco.com**

Internet **www.pasco.com**

Telephone 7:00am - 4:30pm PST

In the U.S. **800.772.8700**

International

1.916.786.3800