
Instruction Sheet for the PASCO Model PS-2525

Adhesive Mylar Foam Covers (10 Pack)

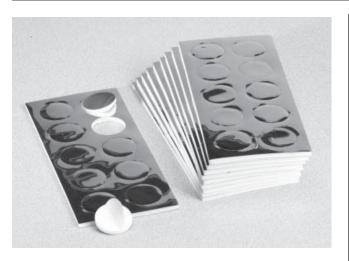

Introduction

The Adhesive Mylar Foam Covers (PS-2525) are replacement units for use with the Fast Response Temperature Probe (PS-2135). The adhesive disks are designed to hold the temperature probe tip and adhere it to a participant's skin or to other flat surfaces. The mylar reduces environmental influences.

Using the Adhesive Mylar Cover

1. Remove one of the adhesive mylar foam cover disks from the sheet set.



2. Carefully remove the protective backing paper.



3. Place the Fast Response Temperature Probe sensing tip in the center of the disk adhesive.



4. Place the adhesive mylar foam disk on a smooth (preferably hairless) skin surface





10101 Foothills Blvd. • Roseville, CA 95747-7100 Phone (916) 786-3800 • FAX (916) 786-8905 • www.pasco.com