
CART ADAPTER ACCESSORY

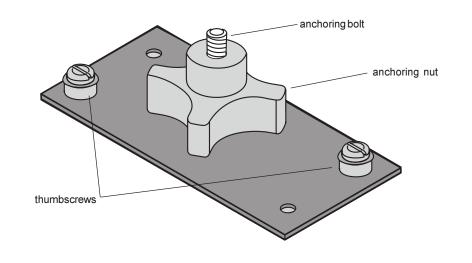

Introduction

The PASCO ME-6743 Cart Adapter Accessory is used to attach the PASCO CI-6742 Motion Sensor II to the PASCO Dynamics Carts and Motorized Carts.

Procedure

- *I*. Attach the plate of the Cart Adapter Accessory with the two thumbscrews (Figure 1).
- 2. Attach the Motion Sensor to the anchoring bolt, turning the Motion Sensor until it is lightly tightened onto the bolt.
- 3. Orient the Motion Sensor in the desired direction.
- *4.* Turn the anchoring nut in a counter-clockwise direction until the Motion Sensor is firmly locked into place.

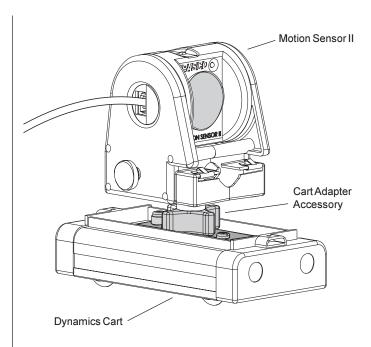


Figure 1

A Motion Sensor attached to a Dynamics Cart with the Cart Adapter Accessory

© 1998 PASCO scientific



10101 Foothills Blvd. • Roseville, CA 95747-7100 Phone (916) 786-3800 • FAX (916) 786-8905 • www.pasco.com



Limited Warranty

PASCO scientific warrants the product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace, at its option, any part of the product which is deemed to be defective in material or workmanship. The warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment after repair will be paid by PASCO scientific.

Address: PASCO scientific 10101 Foothills Blvd. Roseville, CA 95747-7100

Phone:	(916) 786-3800
FAX:	(916) 786-8905
e-mail:	techsupp@pasco.com
web:	www.pasco.com

