
Instruction Manual and Experiment Guide for the PASCO scientific Model CI-6728 012-06619A 10/97

FLUORIDE ION SELECTIVE ELECTRODE

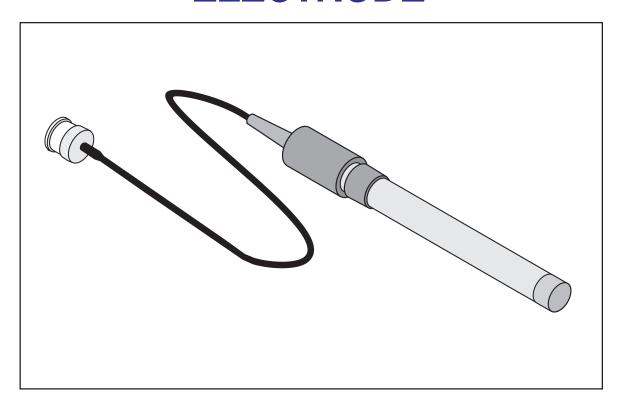

© 1997 PASCO scientific \$7.50







Always use eye protection and gloves when working with chemicals.



Table of Contents

Introduction	1
Theory	1
Equipment	2
Required Solutions	3
General Preparation	4
Electrode Preparation Electrode Slope Check Using <i>Science Workshop</i>	
Measurement	5
Measuring Hints	5
Sample Requirements	
Units of Measurement	5
Measurement Procedure	
Direct Measurement of Fluoride	
Direct Measurement of Fluoride in Water	
Direct Measurement of Fluoride in Acid Solutions	
Direct Measurement of Fluoride in Alkaline Solutions	
Low Level Fluoride Determination	
Titration Procedure for Fluoride Determination	
Electrode Characteristics	
Reproducibility	
Interferences	
Complexation	
Temperature Influences	
Electrode Response	
Limits of Detection	13
pH Effects	13
Electrode Life	14
Maintenance	14
Electrode Storage	14
Polishing the Membrane	14
Specifications	15
Troubleshooting Guide	15
Plastic-ware	
Electrode	15
Standards & Reagents	16
Sample	
Technique	16
Troubleshooting Hints	17
Technical Support	19



Copyright, Warranty, and Equipment Return

Please—Feel free to duplicate this manual subject to the copyright restrictions below.

Copyright Notice

The PASCO scientific 012-06619A manual is copyrighted and all rights reserved. However, permission is granted to non-profit educational institutions for reproduction of any part of the Fluoride Ion Selective Electrode manual providing the reproductions are used only for their laboratories and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Limited Warranty

PASCO scientific warrants the product to be free from defects in materials and workmanship for a period of one year from the date of shipment to the customer. PASCO will repair or replace at its option any part of the product which is deemed to be defective in material or workmanship. The warranty does not cover damage to the product caused by abuse or improper use. Determination of whether a product failure is the result of a manufacturing defect or improper use by the customer shall be made solely by PASCO scientific. Responsibility for the return of equipment for warranty repair belongs to the customer. Equipment must be properly packed to prevent damage and shipped postage or freight prepaid. (Damage caused by improper packing of the equipment for return shipment will not be covered by the warranty.) Shipping costs for returning the equipment after repair will be paid by PASCO scientific.

Credits

Author: Peter Boyle Editor: Steve Miller

Equipment Return

Should the product have to be returned to PASCO scientific for any reason, notify PASCO scientific by letter, phone, or fax BEFORE returning the product. Upon notification, the return authorization and shipping instructions will be promptly issued.

NOTE: NO EQUIPMENT WILL BE ACCEPTED FOR RETURN WITHOUT AN AUTHORIZATION FROM PASCO.

When returning equipment for repair, the units must be packed properly. Carriers will not accept responsibility for damage caused by improper packing. To be certain the unit will not be damaged in shipment, observe the following rules:

- ① The packing carton must be strong enough for the item shipped.
- ② Make certain there are at least two inches of packing material between any point on the apparatus and the inside walls of the carton.
- ③ Make certain that the packing material cannot shift in the box or become compressed, allowing the instrument come in contact with the packing carton.

Address: PASCO scientific

10101 Foothills Blvd. P.O. Box 619011

Roseville, CA 95678-9011

Phone: (916) 786-3800 FAX: (916) 786-3292

email: techsupp@pasco.com web: www.pasco.com

Introduction

The PASCO scientific Fluoride Ion Selective Electrode is used to quickly, simply, accurately, and economically measure fluoride ions in aqueous solutions.

Theory

The PASCO scientific Fluoride Ion Selective Electrode consists of a single crystal of lanthanum fluoride as the membrane, bonded into a glass or an epoxy body. Only fluoride ions are mobile in the ionic conductor crystal. When the membrane comes in contact with a solution containing fluoride ions, a potential develops across the membrane. This electrode potential is measured against a constant reference potential, using an ISE Amplifier and a *ScienceWorkshop* computer interface, and depends on the level of free fluoride ions in the solution. The Nernst equation describes the level of fluoride ions in the solution corresponding to the measured potential:

$$E = E_0 - S \log X$$

where:

E = measured electrode potential

 E_0 = reference potential (a constant)

 $S = \text{electrode slope} \ (\approx 57 \, \frac{\text{mV}}{\text{decade}})$

X = level of fluoride ions in solution

The activity, X, represents the effective concentration of free fluoride ions in the solution. Total fluoride concentration, C_t , may include some bound as well as free fluoride ions. Since the electrode only responds to free ions, the concentration of the free ions, C_t , is found by:

$$C_f = C_t - C_b$$

where C_h represents the concentration of all bound or complexed fluoride ions.

The activity is related to the free ion concentration, C_f , by the activity coefficient, γ , by:

$$X = \gamma C_t$$

Activity coefficients vary, depending on total ionic strength, I, defined as:

$$I = \frac{1}{2} \Sigma C_X Z_{X^2}$$

where:

 C_X = concentration of ion X

 Z_X = charge of ion X

 Σ = sum of all of the types of ions in the solution



In the case of high and constant ionic strength relative to the sensed ion concentration, the activity coefficient, γ , is constant and the activity, X, is directly proportional to the concentration.

All samples and standards containing fluoride ions have TISAB (Total Ionic Strength Adjustment Buffer) added so that the background ionic strength is high and constant relative to variable concentrations of fluoride. The recommended ISA (Ionic Strength Adjuster) for the fluoride electrode is TISAB though similar solutions can be used as long as they do not contain ions that would interfere with the electrode's response to fluoride.

The reference electrode must also be considered. When two solutions of different composition are brought into contact with one another, liquid junction potentials arise. Millivolt potentials occur from the interdiffusion of ions in the two solutions. Electrode charge ,will be carried unequally across the solution boundary resulting in a potential difference between the two solutions, since ions diffuse at different rates. When making measurements, it is important to remember that this potential be the same when the reference is the standardizing solution as well as in the sample solution or the change in liquid junction potential will appear as an error in the measured electrode potential.

The composition of the liquid junction filling solution in the reference electrode is most important. The speed with which the positive and negative ions in the filling solution diffuse into the sample should be equitransferent. No junction potential can result if the rate at which positive and negative charge carried into the sample is equal.

Equipment


Included

- Fluoride Ion Selective Electrode
- Fluoride Ion Selective Electrode fill solution
- pipette for fill solution
- polishing strips

Additional Required:

Required Equipment

- PASCO CI-6738 ISE (Ion Selective Electrode) Amplifier
- ScienceWorkshop 2.2.5 or higher
- PASCO Science Workshop Computer Interface
- Semilogarithmic 4-cycle graph paper for preparing calibration curves when using the meter in the my mode.
- A magnetic stir plate
- Lab-ware made of plastic, not glass

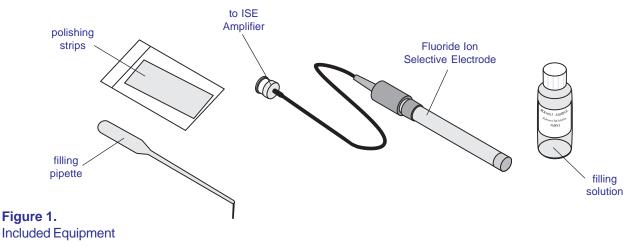

Required Solutions

The stock solutions described in this section may be created as described in the text or ordered directly from PASCO. The solutions available for order, and their respective prices are listed on the 'ISE Working Solution Price List'.

- Deionized or distilled water for solution and standard preparation.
- Fluoride Standard Solution, 0.1 M NaF

To prepare this solution, half fill a one liter volumetric flask with distilled water and add 4.2 grams of reagent-grade sodium fluoride. Swirl the flask gently to dissolve the solid. Fill the flask to the mark with distilled water, cap, and upend several times to mix the solution.

• Fluoride Standard, 1000 ppm F ⁻¹

To prepare this solution, half fill a one liter volumetric flask with distilled water and add 2.21 grams of reagent-grade sodium fluoride. Swirl the flask gently to dissolve the solid. Fill the flask to the mark with distilled water, cap, and upend several times to mix the solution.

• Fluoride Standard, 100 ppm F ⁻¹

To prepare this solution, half fill a one liter volumetric flask with distilled water and add 0.22 grams of reagent-grade sodium fluoride. Swirl the flask gently to dissolve the solid. Fill the flask to the mark with distilled water, cap, and upend several times to mix the solution.

• Total Ionic Strength Adjuster Buffer, TISAB 1

TISAB 1 is used to adjust the pH of the solution, de-complex fluoride and provide a constant background ionic strength. To prepare this solution, half fill a four liter beaker with distilled water, place the beaker on a magnetic stirrer, add a large stirring bar, and begin stirring. Slowly add 230 ml of concentrated acetic acid, 232 grams of reagent-grade sodium chloride, and 16 grams of reagent-grade CDTA. After the solids have dissolved, allow the solution to cool to room temperature. Slowly add 150 grams of reagent-grade sodium hydroxide. After the solids have dissolved, allow the solution to cool to room temperature. Calibrate a pH electrode and adjust the pH to 5.25 with small addition of 5 M NaOH. Fill to the mark with distilled water.

• Low Level Total Ionic Strength Adjuster Buffer, TISAB 2

Use when measuring in samples containing less than 2X10⁻⁵ M (0.4 ppm) fluoride and containing no fluoride complexing agents. To prepare this solution, place about 2000 ml distilled water in a four liter beaker. Add 57 ml glacial acetic acid and 55 grams sodium chloride. Place the beaker on a magnetic stirrer, add a stirring bar and begin stirring. Immerse a calibrated pH electrode into the solution. Slowly add 5 M NaOH until the pH is 5.25. Allow the solution to cool and fill to the mark with distilled water.

• Total Ionic Strength Adjuster Buffer, TISAB 3

TISAB 3 will complex more than 100 ppm of aluminum or iron in the presence of 1 ppm fluoride ion. There will be an error of approximately 5% in the measurement of 1 ppm fluoride in the presence of 200 ppm aluminum or iron. To prepare this solution, add about 2000 ml of distilled water to a four liter beaker. Place the beaker on a magnetic stirrer, add a large stirring magnet, and begin stirring. Slowly add 336 ml of concentrated HCl (36 - 38%), 968 grams of TRIS (hydroxymethylaminomethane) and 920 grams of sodium tartrate ($Na_2C_4H_4O_6$ $2H_2O$). After the solid has dissolved, allow the solution to cool to room temperature. Fill to the mark with distilled water.




General Preparation


Electrode Preparation

- Remove the rubber cap covering the electrode tip and the rubber insert covering the filling hole of the reference electrode. Fill the combination electrode or the reference electrode with the filling solution shipped with the electrode to a level just below the fill hole.
- 2. Connect the Fluoride Ion Selective Electrode to the ISE Amplifier and insert the DIN connector of the ISE Amplifier into analog channel A or B on a PASCO Computer Interface (Figures 2a and 2b).

Electrode Slope Check Using ScienceWorkshop (check electrodes each day)

- 1. To a 150 ml plastic beaker, add 50 ml of distilled water and 50 ml of TISAB. Place the beaker on a magnetic stirrer and begin stirring at a constant rate. Start the *ScienceWorkshop* software, select the Ion Selective Electrode sensor, open a Digital display, change the number of digits to the right of the decimal from 1 to 3, and begin monitoring data. Lower the electrode tip into the solution.
- 2. Using a pipette, add 1 ml of 0.1 M, 1000 ppm, or 100 ppm fluoride standard to the beaker. When the reading has stabilized, record the Voltage reading in the Digits display.
- 3. Using a pipette, add 10 ml of the same fluoride standard used above to the beaker. When the reading has stabilized, record the Voltage reading in the Digits display.
- 4. Determine the difference between the two readings. The electrode is operating correctly if the potential has changed by 57±2mV, assuming the solution temperature is between 20 °C and 25 °C. See the *Troubleshooting* section if the potential change is not within this range.
 - Note: Slope is defined as the change in potential observed when the concentration changes by a factor of 10.

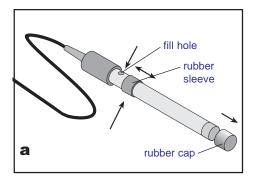

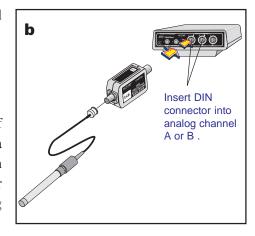

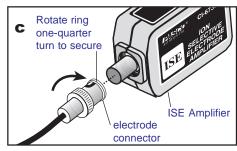


Figure 2
Equipment Setup. a: filling the electrode with filling solution; b & c: connecting the electrode to the ISE Amplifier and to the computer interface



Measurement

Measuring Hints

• All samples and standards should be at the same temperature for precise measurement. A difference of 1 °C in temperature will result in a 2% measurement error.

- Constant, but not violent, stirring is necessary for accurate measurement. Magnetic stirrers can generate sufficient heat to change the solution temperature. To counteract this effect, place a piece of insulating material, such as a styrofoam sheet, between the stirrer and the beaker.
- Always rinse the electrodes with distilled water and blot dry between measurements. Use a clean, dry tissue to prevent cross-contamination.
- For samples with high ionic strength, prepare standards whose composition is similar to the sample.
- Always check to see that the membrane is free from air bubbles after immersion into standard or sample.

Sample Requirements

- All samples must be aqueous and not contain organics which can dissolve the epoxy electrode body and/
 or the cement bonding the sensing crystal to the electrode body. Inorganic solutions will not effect the
 electrode. Infrequent measurements in solutions containing methanol, acetone, or dioxane are permitted.
 Highly polar solvents, such as CHC1₃ or DMF, should not be contained in the samples.
- The addition of TISAB to samples and standards will adjust the pH to 5.0 5.5. Samples must be above pH 5 to avoid forming complexes with hydrogen ions and below pH 7 to avoid interference by hydroxide ions.
- The temperature of the standard solutions and of the sample solutions should be the same and below $80\,^{\circ}\text{C}$.
- The use of TISAB 1 also preferentially forms complexes with aluminum and with iron, breaking the complexes that fluoride forms with these ions. With 1 ppm fluoride present, up to 3 5 ppm aluminum or iron is complexed. If higher levels of aluminum or iron are present, use TISAB 3.

Units of Measurement

Fluoride concentrations are measured in units of ppm as fluoride, moles per liter, or any other convenient concentration unit. Table 1 indicates some concentration units and conversion factors.

TABLE 1: Concentration Unit Conversion Factors

ppm F ⁻¹	moles/liter
190.00	$1.0 \times 10^{-2} M$
19.00	$1.0 \times 10^{-3} M$
1.90	$1.0 \times 10^{-4} M$



Measurement Procedure

Direct Measurement

Direct measurement is a simple procedure for measuring a large number of samples. A single reading is all that is required for each sample. The ionic strength of samples and standards should be made the same by adjustment with TISAB for all fluoride solutions. The temperature of both sample solution and standard solution should be made the same.

Direct Measurement of Fluoride

Note: A calibration curve is constructed on semilogarithmic paper. The measured electrode potential (linear axis) is plotted against the standard concentration (log axis). In the <u>linear region of the curve</u>, only two standards are necessary to determine a calibration curve. Calibration solutions close to the anticipated value of the "unknown" should be chosen. In the <u>nonlinear region</u>, additional points must be measured. The direct measurement procedures given are for the linear portion of the curve. The nonlinear portion of the curve requires the use of low level procedures.

- 1. By serial dilution, prepare three standard solutions from the 0.1 M, 1000 ppm, or the 100 ppm stock standard. The resultant concentrations should be 10⁻² M, 10⁻³ M, and 10⁻⁴ M or 100, 10, and 1 ppm. Add 50 ml of TISAB 1 or TISAB 3 to each 50 ml of standard. When Calibrating, assume that the added TISAB has no effect on the standard concentration.
- Place the most dilute solution on the magnetic stirrer and begin stirring at a constant rate. After assuring
 that *ScienceWorkshop* is operating, lower the electrode tip into the solution. When the reading has
 stabilized, record the voltage reading indicated in the Digits display.
- 3. Place the mid-range solution on the magnetic stirrer and begin stirring. After rinsing the electrodes with distilled water, blot dry, and immerse the electrodes in the solution. When the reading has stabilized, record the voltage reading indicated in the Digits display.
- 4. Place the most concentrated solution on the magnetic stirrer and begin stirring. After rinsing the electrodes with distilled water, blot dry, and immerse the electrodes in the solution. When the reading has stabilized, record the voltage reading indicated in the Digits display.
- 5. Using the semilogarithmic graph paper, plot the voltage reading (linear axis) against the concentration. (log axis). Extrapolate the curve down to about 1.0Xl0 ⁻⁵ M. A typical calibration curve can be found in Figure 3.



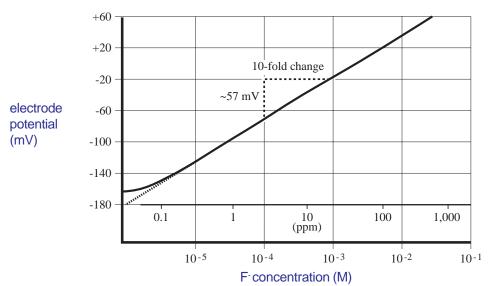


Figure 3Typical Fluoride Ion Selective Electrode calibration curve

- 6. To a clean, dry, 150 ml plastic beaker, add 50 ml of sample and 50 ml of TISAB 1 or TISAB 3. Place the beaker on the magnetic stirrer and begin stirring. Rinse the electrodes with distilled water, blot dry, and lower the electrode tip into the solution. When the reading has stabilized, record the Voltage reading in the Digits display. Using the calibration curve, determine the sample concentration.
- 7. The calibration should be checked every 1 2 hours. Assuming no change in ambient temperature, place the electrode tip in the mid-range standard. After the reading has stabilized, compare it to the original reading recorded in Step 3 above. A reading differing by more than 0.5 mV or a change in the ambient temperature will necessitate the repetition of Steps 2 5 above. A new calibration curve should be prepared daily.

Direct Measurement of Fluoride in Water (ASTM D1179, Method B)

The procedure is used to determine total fluoride concentration in water in ppm. TISAB is added to standards and to samples in order to break fluoride complexes of iron and aluminum, adjust the pH, and provide a constant ionic strength. Fluoride sample concentration is determined, using this method, independent of the level or nature of dissolved minerals.

- 1. Prepare 2, 1, and 0.5 ppm fluoride standards by serial dilution of the 100 ppm fluoride standard.
- 2. To each 50 ml of standard, add 50 ml of TISAB 1 or TISAB 3.
- 3. Calibrate the meter as previously described by using the 2, 1, and 0.5 ppm standards. The calibration curve should be drawn on semi-logarithmic 2-cycle graph paper or *ScienceWorkshop*.

Direct Measurement of Fluoride in Acid Solutions

Hydrogen ion complexes a portion of the fluoride ion in solutions with a pH below 5, forming HF or HF₂⁻¹, which cannot be detected by the electrodes. Adjustment to weakly acidic/weakly basic range before making the fluoride determination is necessary, but not with strongly basic solutions. The use of sodium acetate buffers the pH above 5 and helps fix the total ionic strength of standards and samples to the same level.



 Dissolve reagent grade sodium acetate (CH₃COONa) in distilled water to prepare a 15% solution. Prepare a sufficient quantity to dilute all standards and samples.

- 2. Prepare a blank solution containing all components of the sample except fluoride. This solution will be used to prepare standards.
- 3. Add fluoride to the blank solution to prepare standards in the concentration range of the unknown solutions. Prepare three standards and a calibration curve as previously described. Add 9 parts of sodium acetate to each 1 part of standard. Fresh standards should be prepared every two weeks if the standard contains less than 10 ppm fluoride.
- 4. Calibrate the electrodes as described in the section Electrode Slope Check.
- 5. After diluting each unknown sample 10:1 with sodium acetate (as in step 3 above), measure the electrode potential and determine the fluoride concentration.

Direct Measurement of Fluoride in Alkaline Solutions

Hydroxide ions interfere with fluoride measurements in basic solutions with a low fluoride content. At a fluoride concentration less than 1.0×10^{-4} M and at a pH of 9.5 or above, the electrode potential reading, caused by the concentration of both hydroxide and fluoride ions, is higher than it would be if fluoride alone were present. (See section entitled *pH Effects*.)

Using a 4 M buffered potassium acetate solution to adjust the pH to between 5 and 6 eliminates hydroxide ion error and raises the total ionic strength of both standards and samples to the same value. The fluoride ion concentration can be determined in the usual manner after both standards and samples are diluted 10:1 with the buffer solution.

- 1. To prepare a 4 M buffered potassium acetate solution, add one part of distilled water, slowly, to two parts of 6 M acetic acid, CH₃COOH, in a large beaker surrounded by a water bath. Slowly add 50% KOH solution to the acetic acid mixture, with constant stirring, until a pH of 5 is reached. Prepare enough buffer to dilute all standards and samples 10:1.
- 2. Following the directions given in the preceding section (*Determination of Fluoride in Acid Solutions*), prepare standards, calibrate the electrodes and measure the unknown samples.

Low Level Fluoride Determination

Use the following low level fluoride measurement procedure in the non-linear portion of the calibration curve. (See Figure 1.) This procedure is used for fluoride samples containing less than $2X10^{-5}$ M or 0.4 ppm fluoride and containing no fluoride complexing agents. Use low level TISAB 2 for both samples and standards. A longer response time should be expected for low level fluoride measurements.

- 1. By serial dilution, prepare a 1.0X10⁻³ M or 10 ppm fluoride standard by diluting the 0.1 M, 1000 ppm, or 100 ppm standard solution. Add 50 ml low level TISAB 2 to 50 ml of standard solution.
- 2. Using a 150 ml beaker, add 50 ml of distilled water and 50 ml of low level TISAB 2. Place the beaker on the magnetic stirrer and begin stirring at a constant rate. Lower the electrode tip into the solution. Start the *ScienceWorkshop* software, select the Ion Selective Electrode sensor, open a Digital display, and begin monitoring data.



3. Increments of the standard should be added to the beaker according to the steps outlined in Table 2 below. After the reading stabilizes, record the Voltage reading in the Digits display after each addition.

TABLE 2: Low Level Measurement Calibration Curve

	Added		Concentration	
Step	Pipette	Volume (ml)	ppm	M
1	A	0.1	0.01	1.0 X 10 ⁻⁶
2	A	0.1	0.02	2.0 X 10 ⁻⁶
3	A	0.2	0.04	4.0 X 10 ⁻⁶
4	A	0.2	0.06	6.0 X 10 ⁻⁶
5	A	0.4	0.10	1.0 X 10 ⁻⁵
6	В	2.0	0.29	2.9 X 10 ⁻⁵
7	В	2.0	0.48	4.8 X 10 ⁻⁵

Pipette A = 1 ml graduated pipette

Pipette B = 2 ml pipette

Solutions: additions of standard/TISAB 2 to 50 ml of distilled water and 50 ml of low level TISAB 2.

- 4. On semi-logarithmic graph paper, plot the concentration (log axis) against the millivolt reading (linear axis) as in Figure 1. Keep the final solution for checking the electrode calibration.
- 5. To a 150 ml plastic beaker, add 50 ml of sample and 50 ml of low level TISAB 2. Place the beaker on a magnetic stirrer and begin stirring. After rinsing the electrode and blotting dry, place the electrode tip into the solution. After stabilization of the reading, read the electrode potential and determine the concentration from the calibration curve. A new low level calibration curve should be prepared daily using fresh standards.

Titration

Titration is a very accurate determination of fluoride. The PASCO Fluoride Ion Selective Electrode can be used as a highly sensitive endpoint detector for titrations of a fluoride-containing sample. Though titrations are more time consuming than direct ion measurements, the results are more accurate and reproducible. Titrations accurate to $\pm 0.2\%$ of the total fluoride concentration of the sample can be performed using lanthanum nitrate as the titrant. Total fluoride concentration should be at least 1.0×10^{-3} M for endpoint detection. Low results are given if aluminum, iron, or trivalent chromium are present at a level of 1% or higher.

Special titration procedures for aluminum, lithium, lanthanum, and thorium also makes use of the fluoride electrode as an endpoint indicator.

Titration Procedure for Fluoride Determination

Dissolve 43.3 grams of reagent grade lanthanum nitrate, La(NO₃)₃• 6H₂O, in about 500 ml distilled water in a 1 liter volumetric plastic flask. Fill to the mark with distilled water. This 0.10 M lanthanum solution will be used for all titrations.



2. Using the 0.1 M fluoride standard, standardize the lanthanum nitrate by titration. To a 150 ml plastic beaker, add approximately 9.0 ml of fluoride standard (accurately measured) and about 50 ml of distilled water. Place the beaker on the magnetic stirrer and begin stirring. Lower the electrode tip into the solution.

- 3. Using a 10 ml plastic burette, add the La (NO₃)₃ titrant in 0.5 1.0 ml increments. Record the Voltage reading in the Digits display against the volume of titrant added. As the electrode potential change increases, add smaller increments, down to 0.1 0.2 ml increments. Continue to add titrant and record the electrode potential against the volume until little change is noted in the Voltage reading even when adding 0.5 1.0 ml increments.
- 4. Using linear graph paper, plot the Voltage reading (y-axis) against the volume (x-axis). The end-point is determined as the steepest slope on the titration curve. Record the endpoint.

 V^0_{t}

- 5. To a 150 ml plastic beaker, add approximately 9.0 ml of sample solution (accurately measured) and about 50 ml of distilled water. Place the beaker on a magnetic stirrer and begin stirring. Lower the rinsed, dried electrode tip in the solution.
- 6. Titrate the sample as in step 3 above. The endpoint is denoted as:

 V^{X}

7. Calculate the sample concentration, C^{x}

$$C_{S}^{X} = \frac{V_{t}^{X} V_{f}^{0}}{V_{t}^{0} V_{f}^{X}} - 2C_{S}^{0}$$

where:

 $C_{s}^{X} = concentration of sample$

 $C_{S}^{0} = fluoride standard concentration$

 V_{t}^{X} = volume of titrant added to achieve the endpoint of unknown sample


 V_{t}^{0} = volume of titrant added to achieve the endpoint in standardization

 V_{f}^{X} = volume of sample used in sample titration

 V_{f}^{0} = volume of standard used in standardization titration



A typical titration curve is shown in Figure 4.

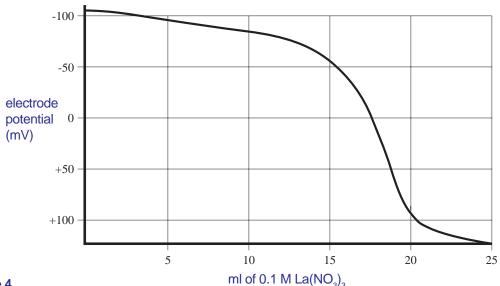

Titration of 25 ml of 0.112 M F $^{-}$ with 0.100 M La (NO $_{3}$) $_{3}$

Electrode Characteristics

Reproducibility

Electrode measurements reproducible to $\pm 2\%$ can be obtained if the electrode is calibrated every hour. Factors such as temperature fluctuations, drift, noise, and variations in illumination limit reproducibility. Reproducibility is independent of concentration within the electrode's operating range.

Interferences

The hydroxide ion, OH $^{-1}$, is an electrode interference. Anions which make the sample more basic, such as ${\rm CO_3}^{-2}$ or ${\rm PO_4}^{-3}$ would increase the OH $^{-1}$ interference, but do not interfere with direct electrode operation. Other anions commonly associated with fluoride, such as Cl $^{-1}$, Br $^{-1}$, I $^{-1}$, SO $_4^{-2}$, HCO $_3^{-1}$, NO $_3^{-1}$ and acetate, do not interfere with correct electrode operation. Most cations do not interfere with the response of the fluoride electrode to fluoride ion.

Complexation

Hydrogen ion, as well as some other multivalent cations, aluminum, silicon, iron+3, will form complexes with fluoride. The total ionic strength of the solution, the pH of the solution, the total fluoride concentration, and the concentration of the complexing agent all contribute to the degree of complexation. TISAB 1 and TISAB 2 complex about 5 ppm aluminum or iron in a 1 ppm fluoride solution. TISAB 3 complexes higher levels of iron and aluminum.



Temperature Influences

Samples and standards should be within $\pm 1\,^{\circ}$ C of each other, since electrode potentials are influenced by changes in temperature. Because of solubility equilibria on which the electrode depends, the absolute potential of the reference electrode changes slowly with temperature. The slope of the electrode, as indicated by the factor "S" in the Nernst equation, also varies with temperature. Table 3 gives values for the 'S" factor in the Nernst equation for the fluoride ion.

TABLE 3: Temperature vs. Theoretical Values for the Electrode Slope

Temperature (°C)	"S"
0	54.20
10	56.18
20	58.16
25	59.16
30	60.15
40	62.13
50	64.11

The temperature range for the PASCO scientific Fluoride Ion Selective Electrode is 0 °C - 80 °C, provided that temperature equilibrium has occurred. Only intermittent use is recommended at temperatures from 80 °C - 100 °C. If the temperature varies substantially from room temperature, equilibrium times up to one hour are recommended.

Electrode Response

Plotting the electrode potential against the fluoride concentration on semi-logarithmic paper results in a straight line with a slope of about 57 mV per decade. (Refer to Figure 1.)

The time needed to reach 99% of the stable electrode potential reading, the electrode response time, varies from one minute or less in highly concentrated solutions to several minutes near the detection limit. See Figure 5.

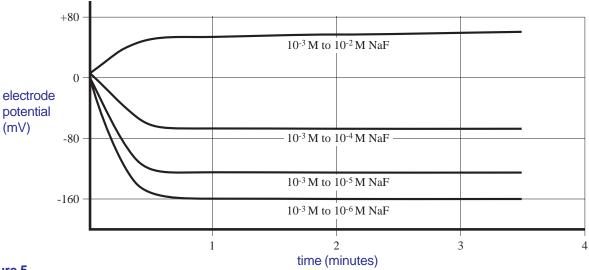


Figure 5Typical electrode time response to step changes in NaF

A drifting potential reading or a decrease in electrode slope may mean that the electrode membrane needs polishing.

Limits of Detection

Fluoride concentration down to $1.0 \times 10^{-6} \,\mathrm{M}$ (0.02 ppm) fluoride can be measured in neutral solutions. Since sample contamination can be a factor in low level fluoride measurements, care must be taken in making determinations below $1.0 \times 10^{-5} \,\mathrm{M}$. The upper limit of detection is a saturated fluoride solution.

pH Effects

Hydrogen complexes a portion of fluoride in solution, forming the un-dissociated acid HF and the ion HF_2^{-1} in acid solutions with a pH below 5. The proportion of free fluoride ion in acid solutions is shown in Figure 6.

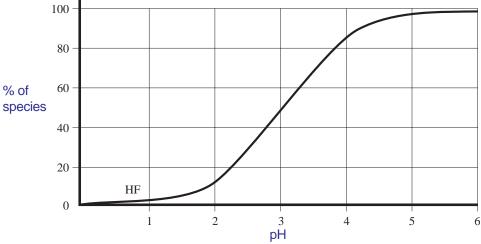


Figure 6Fraction of free F - as a function of solution pH, where the hydrogen is the only complexing agent

When the level of hydroxide is greater than one-tenth the level of fluoride ion present, hydroxide ion interferes with electrode response to fluoride. As an example, no hydroxide interference with fluoride measurements take place at pH 7 when the hydroxide concentration is 1.0×10^{-7} or less. As the pH increases, the hydroxide interference becomes appreciable. At pH 10, the hydroxide ion concentration is 1.0×10^{-4} M and no error is found in measurements of 1.0×10^{-2} M fluoride. At the same hydroxide ion concentration and a fluoride concentration of 1.0×10^{-4} M, about a 10% measurement error appears. At a fluoride concentration of 1.0×10^{-5} M, considerable error exists in a pH 10 solution. Figure 7 illustrates these errors.

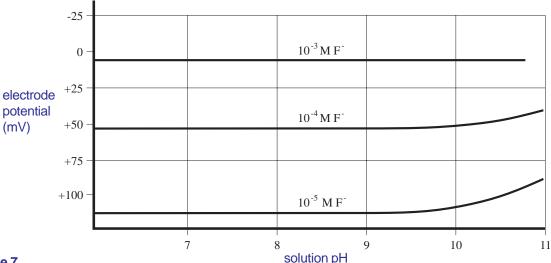


Figure 7 Electrode response in alkaline solution



The addition of TISAB 1 or VISAS 2 to all fluoride samples and standards buffers the pH between 5.0 - 5.5 to help avoid hydroxide interferences or the formation of hydrogen complexes of fluoride. VISAB 3 adjusts the pH to about 8.5 and should not be used for very low level measurements.

Electrode Life

The fluoride electrode will last six months in normal laboratory use. On-line measurements might shorten operational lifetime to several months. In time, the response time will increase and the calibration slope will decrease to the point calibration is difficult and electrode replacement is required.

Maintenance

Electrode Storage

The fluoride electrode may be stored for short periods of time in 1.0×10^{-2} M fluoride solution with TISAB added. For longer storage (longer than two weeks), rinse and dry the sensing pellet and cover the membrane tip with any protective cap shipped with the electrode. The reference portion of the combination electrode should be drained of filling solution, if refillable, and the rubber insert placed over the filling hole. The fluoride electrode should never be stored in distilled water.

Polishing the Membrane

- 1. If using polishing paper, cut off a 1 2" piece and place it face up on the lab bench.
- 2. Put a few drops of distilled or deionized water in the center of the paper.
- 3. Holding the paper (cotton) steady with one hand, bring the membrane of the electrode down perpendicular to the paper and, with a slight swirling motion, gently polish the tip of the electrode against the surface of the polishing paper (cotton) for a few seconds.
- 4. Rinse the electrode surface with distilled or deionized water and soak the electrode tip in standard solution for about five minutes before use.
- 5. If using jeweller's rouge, place a cotton ball on the table top and flatten it using the bottom of a beaker.
- 6. Put 1 2 drops of distilled or deionized water in the center of the cotton pad.
- 7. Add a small amount of jeweller's rouge to the damp cotton.
- 8. Continue with Steps 3 and 4 above.



Specifications

Concentration Range: saturated solutions to 1.0X10⁻⁶ M (0.02 ppm)

pH Range: 5 to 7 at 1.0×10^{-6} M F $^{-1}$ (0.02 ppm F $^{-1}$)

5 to 11 at $1.0 \times 10^{-1} \,\mathrm{M}\,\mathrm{F}^{-1} (1900 \,\mathrm{ppm}\,\mathrm{F}^{-1})$

Temperature Range: $0 \,^{\circ}\text{C} - 80 \,^{\circ}\text{C}$

Resistance: 150 to 200 kohm

Reproducibility: $\pm 2\%$

Storage: in fluoride standard with TISAB added

Size: length = 110 mm

diameter = 12 mmcable length = 1 m

Troubleshooting Guide

The goal of troubleshooting is the isolation of a problem through checking each of the system components in turn: the meter, the plastic-ware, the electrodes, the standards & reagents, the sample, and the technique.

Plastic-ware

Clean plastic-ware is essential for good measurement. Be sure to wash the plastic-ware well with a mild detergent and rinse very well with distilled or deionized water.

Electrode

The electrode may be checked by using the procedure found in the sections entitled *Electrode Slope Check*.

- 1. Be sure to use distilled or deionized water when following the procedures given in *Electrode Slope Check*.
- 2. If the electrode fails to respond as expected, see the sections *Measuring Hints* and *Electrode Response*. Repeat the slope check.
- 3. If the electrodes still fail to respond as expected, substitute another Fluoride Ion Selective Electrode (if available) that is known to be in good working order for the questionable electrode.
- 4. If the problem persists, the reagent may be of poor quality, interferences in the sample may be present or the technique may be faulty. (See *Standards & Reagents*, and *Technique* sections below.)



5. If another electrode is not available for test purposes, or if the electrode in use is suspect, review the instruction manual and be sure to:

- Clean and rinse the electrodes thoroughly.
- Prepare the electrodes properly.
- Use the proper filling solution.
- Adjust the pH and the ionic strength of the solution by the use of the proper TISAB.
- Measure correctly and accurately.
- Review Troubleshooting Hints.

Standards & Reagents

Whenever problems arise with the measuring procedure that has been used successfully in the past, be sure to check the standard and reagent solutions. If in doubt about the credibility of any of the solutions, prepare them again. Errors may result from contamination of the TISAB, incorrect dilution of standards, poor quality distilled/deionized water, or a simple mathematical miscalculation.

Sample

Look for possible interferences, complexing agents, or substances which could affect the response or physically damage the sensing electrode if the electrode works perfectly in the standard, but not in the sample.

Try to determine the composition of the samples prior to testing to eliminate a problem before it starts. (See *Measuring Hints*, *Sample Requirements*, and *Interferences*.)

Technique

Be sure that the electrode's limit of detection has not been exceeded. Be sure that the analysis method is clearly understood and incompatible with the sample.

Refer to the instruction manual again. Reread General Preparation and Electrode Characteristics.

If trouble still persists, call PASCO scientific Technical Support.



Troubleshooting Hints

Symptom	Possible Causes	Next Step	
Out of Range Reading	defective electrode	check electrode operation	
	electrodes not plugged in properly	unplug electrodes and reseat electrodes	
	reference electrode not filled	replenish reference filling solution	
	electrodes not in solution	put electrodes in solution	
	air bubble on membrane	remove bubble by re-dipping electrode	
Noisy or Unstable Readings (readings continuously or rapidly	electrode exposed to interferences	soak electrode in fluoride standard with added TISAB	
changing)	defective electrode	replace electrode	
	TISAB not used	use recommended TISAB	
	stirrer not grounded	ground stirrer	
	air bubble on membrane	remove bubble by re-dipping electrode	
Drift (reading slowly changing in one direction)	samples and standards at different temperatures	allow solutions to come to room temperature before measurement	
	electrode exposed to complexing agents	check section entitled Complexation	
	incorrect reference filling solution	use recommended filling solution	



Symptom	Possible Causes	Next Step	
Low Slope or No Slope	standards contaminated or incorrectly made	prepare fresh standards	
	TISAB not used	use recommended TISAB	
	standard used as TISAB	use TISAB	
	electrode exposed to complexing agents	check section entitled Complexation	
	air bubble on membrane	remove bubble by re-dipping probe	
"Incorrect Answer" (but calibration curve is good)	incorrect scaling of semilog paper	plot voltage potential on the linear axis. On the log axis, be sure concentration numbers within each decade are increasing with increasing concentration.	
	incorrect sign	be sure to note sign of millivolt reading correctly	
	incorrect standards	prepare fresh standards	
	wrong units used	apply correct conversion factor: $1.0 \times 10^{-3} \mathrm{M} = 19.0 \mathrm{ppm} \;\mathrm{as}\;\mathrm{F}^{-1}$	
	electrode exposed to complexing agents	check section entitled <i>Complexation</i>	



Technical Support

Feedback

If you have any comments about the product or manual, please let us know. If you have any suggestions on alternate experiments or find a problem in the manual, please tell us. PASCO appreciates any customer feedback. Your input helps us evaluate and improve our product.

To Reach PASCO

For technical support, call us at 1-800-772-8700 (toll-free within the U.S.) or (916) 786-3800.

fax: (916) 786-3292

e-mail: techsupp@pasco.com

web: www.pasco.com

Contacting Technical Support

Before you call the PASCO Technical Support staff, it would be helpful to prepare the following information:

- ➤ If your problem is with the PASCO apparatus, note:
 - Title and model number (usually listed on the label);
 - Approximate age of apparatus;
 - A detailed description of the problem/ sequence of events (in case you can't call PASCO right away, you won't lose valuable data);
 - If possible, have the apparatus within reach when calling to facilitate description of individual parts.
- ➤ If your problem relates to the instruction manual, note:
 - Part number and revision (listed by month and year on the front cover);
 - Have the manual at hand to discuss your questions.