


# Force Platform Handle Set

PS-2548

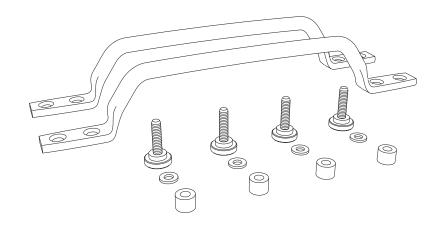

Additional Equipment Required	Part Number
Force Platform (1D)	PS-2141

### Introduction

The PASCO model PS-2548 Force Platform Handle Set is designed as an accessory to the PASCO Force Platform. The handles may be attached to the top of the Force Platform, the bottom of the Force Platform, or both.

# **Experiments**


Please refer to the experiment guide that comes with the applicable PASCO equipment for specific experiments.

### Installation procedure

#### Installation on the top of the Force Platform

1. The corners of the Force Platform are identified by numbers molded into the top of the case. Place one handle spanning corners 1-4. Place the other handle spanning corners 2-3.


2. Use the supplied thumb screws and washers to firmly attach the handles as shown in the following illustration. The metal spacers supplied with the thumb screws are used when attaching handles to the bottom of the Force Platform and should not be used when installing handles on the top.

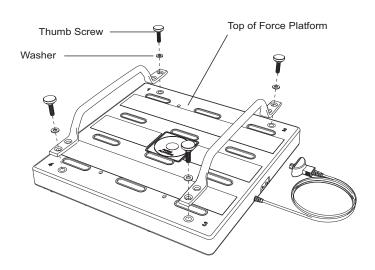

Handle installation on top of Force Platform

#### Installation on the bottom of the Force Platform

- 1. With the bottom of the Force Platform facing up, unscrew the four feet and set them aside.
- 2. Place a spacer in the recess at each attachment point.
- 3. Use the thumb screws with washers to firmly attach the handles to the Force Platform as shown in the following illustration.

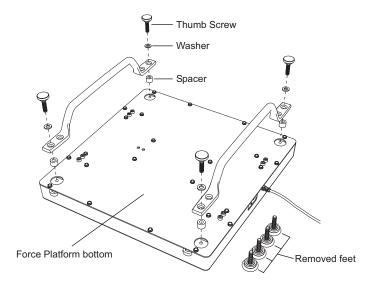

Handle installation on bottom of Force Platform

## **Technical Support**

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific

10101 Foothills Blvd. Roseville, CA 95747-7100

Phone: 916-786-3800 (worldwide)

800-772-8700 (U.S)

Fax: (916) 786-3292
Web: www.pasco.com
Email: support@pasco.com

#### **Limited Warranty**

For a description of the product warranty, see the PASCO catalog.

#### Copyright

The PASCO scientific 012-09470B Force Platform Handle Set Instruction Sheet is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

#### Trademarks

PASCO, PASCO scientific, DataStudio, PASPORT, and ScienceWorkshop are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.