




PASCO Structures Systems Large Structures Set ME-7003

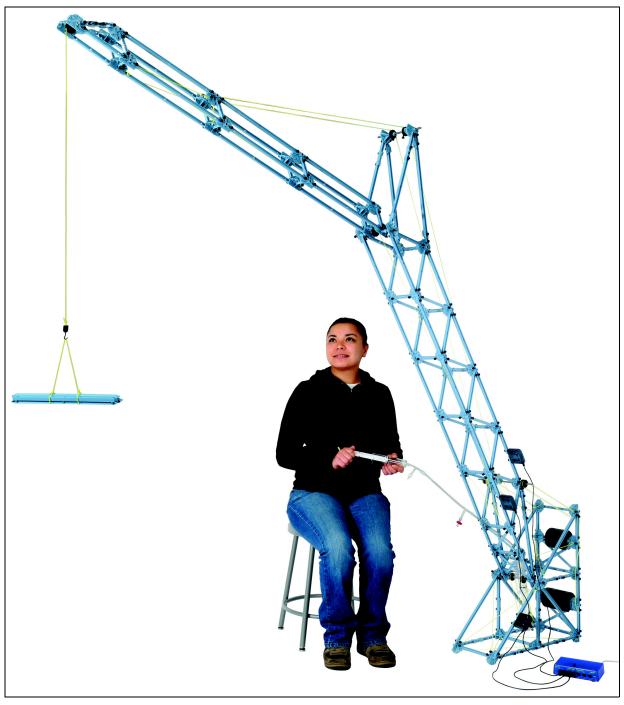

800-772-8700

www.pasco.com

The cover page shows an Angle Crane model built with components from the ME-7003 Large Structures Set. The picture shows Load Cells, a Load Cell Amplifier, Large Slotted Masses, and the Hydraulics Structures Set which are available separately. The ME-7003 Large Structures Set can be used to build over a dozen different models. Instructions for building twenty models are included in this manual.

Table of Contents

Included Items
Recommended Equipment
Introduction
About the Components
Operation
Adding Load Cells
Large Structures Construction
1. Car Building Jump
2. Cable Stayed Bridge
3. Arch Truss Bridge
4. Roller Coaster
5. Canyon Car Jump
6. Suspension Bridge
7. High Road/Low Road
8. Tied Arch Bridge
9. Baltimore Bridge
10. Double Tied Arch Bridge
11. Cantilever Truss Bridge
12. Brachistochrone
13. Drawbridge
14. Vertical Lift Drawbridge
15. Tower Crane
16. Arch Causeway
17. Skyscraper
18. House Frame
19. Angle Crane

20. Windmill	50
Technical Support, Warranty, and Copyright	54
Windmill Sail Template	55

Large Structures Set

ME-7003



Included Items	Qty	Included Items	Qty
ME-6974 Large Structures Parts	1 set	ME-6998A Axle Spares	1 set
ME-6986 Structures Rod Clamps (2/set)	1 set	ME-6999A Angle Connectors	1 set
ME-6985 Flexible I-Beams	1 set	ME-7002 Connector Spares	2 sets
ME-6987 Flat Structures Members	1 set	ME-7008 #6 I-Beam Spares	1 set
ME-6988A Force Platform Structure	1 set	ME-9814 Coaster Track (9.1 m)	1
ME-6993 Truss Set Members	3 sets	ME-9839 Mini-car, Green	1
ME-6994 Truss Set Screws (75/set)	6 sets	ME-9840 Mini-car, Yellow	1
ME-6996 Cord Lock Spares	1 set	ME-9856 Mini-car Starter Bracket	1
ME-6997 Full Round Connector Set	1 set	Container (not shown)	2

The ME-7003 Large Structures Set can be used with other sets from the PASCO Structures System* such as the following:.

ME-6984 Hydraulics Structures Set	ME-6991 Bridge Set
ME-6990 Truss Set	ME-6992A Advanced Structures Set

*See the PASCO catalog or Web site at www.pasco.com for information about the PASCO Structures System.

The following equipment* is recommended for use with a PASCO Interface to measure tension and compression forces in the structures or to measure the motion of a Mini Car as it moves on a structure.

Recommended Equipment*	Recommended Equipment*		
Load Cell Amplifier (PS-2198, PS-2206 or CI-6464)	PS-2204 Displacement Sensor		
PS-2200 100 N Load Cell	PS-2159 PASPORT Digital Adapter		
PS-2201 5 N Load Cell	ME-9498A Photogate Head		

*See the PASCO catalog or Web site at www.pasco.com for more information about the recommended items.

Introduction

The ME-7003 Large Structures Set is one part of the PASCO Structures System and can be combined with other parts of the Structures System. The Large Structures Set allows you to build a variety of realistic truss structures. You can add Load Cells to the structures in order to measure the forces anywhere in the structure. This manual describes the construction of sixteen structures.

Load Cell Amplifier (PS-2198 or CI-6464) - The amplifier requires a PASPORT or *ScienceWorkshop* interface to connect to a computer. The PS-2198 accepts up to six Load Cells and the CI-6464 supports one Load Cell.

100 N Load Cell (PS-2200) and 5 N Load Cell (2201) - Strain gauges mounted on a beam with no electronics so a Load Cell requires a Load Cell Amplifier (PS-2198), Dual Load Cell Amplifier (PS-2206), *ScienceWorkshop* Load Cell Amplifier (CI-6464), or CI Sensor Voltage Monitor (CI-6611).


Displacement Sensor (PS-2205) - A PASPORT Sensor and a digital displacement indicator designed to measure the deflection of parts of a structure such as a truss or a bridge as forces are applied.



About the Components

Large Structures Parts (ME-6974)

The ME-6974 Large Structures Parts set consists of two Mini Car Photogate Flags, two Couplers for the Coaster Track, and a set of 24 Track Clips. The Photogate Flags mount on any Mini Car. When the flag on a moving car travels through a photogate and interrupts the beam, the car's motion can be measured. The Couplers are used to connect Coaster Tracks together, and the Track Clips are used to connect a Coaster Track to a PASCO structure.)

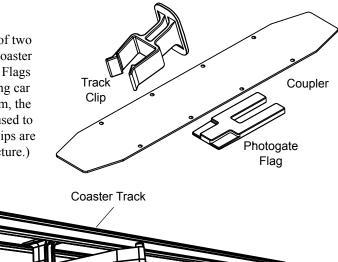

Beam



Channel

Rotate the clip one-quarter turn.

Use track clips to connect the Coaster Track to the structure beams. Slide the clip into the channel on the underside of the track. Move the clip into position and lock it into place with a quarter twist as shown.

Flexible I-Beams (ME-6985)


The Flexible I-Beam Set contains forty-six flexible beams of three different lengths. The flexible beams have the same lengths as three of the I-beams from the Truss Set Members (ME-6993). The flexible beams allow you to explore the concept of structural failure. By replacing a #3, #4, or #5 I-beam in a structure with one of the flexible I-beams you can observe the effect of a load that causes deformation. These I-beams are designed to demonstrate stress and failure and then return to their original shape once the load is removed.

Flexible I-Beam Set (ME-6985)	Qty	Equivalent Truss Set Members
#3 Flexible I-Beam (11.5 cm)	18	#3 I-Beam
#4 Flexible I-Beam (17 cm)	18	#4 I-Beam
#5 Flexible I-Beam (24 cm)	10	#5 I-Beam

Please note that the flexible I-beams are a different color (beige) than the other I-beams (light blue).


Structures Rod Clamp (ME-6986)

Use a rod clamp to attach a structure beam to any 12.7 mm (1/2") diameter rod.

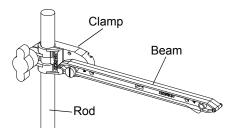

Flat Structures Members (ME-6987)

This set contains three types of flat structures: Flat 3 X 4 Beam (19 cm), Flat #4 Beam (17 cm), and Flat 2 X 3 Beam (12.5 cm). There are sixteen of each type of beam in the set.

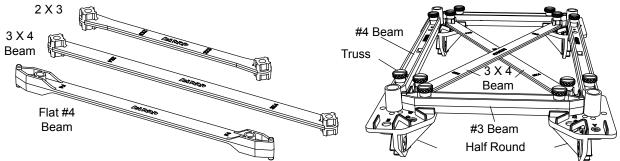

Force Platform Structure (ME-6988A)

The PASCO model ME-7003 Force Platform Structures Bracket includes two brackets and four thumbscrews. The adapter bracket is designed to connect members of the PASCO Structures System to a PASCO Force Platform (not included). The brackets can also serve as foundation plates for larger structure models. (Please see the Force Platform Structures Bracket instruction sheet for more information.)

Truss Set Members (ME-6993)


The Truss Set Members consists of five types of "I-beams" and one type of connector. Each set has the following quantities:

Item	Qty	Item	Qty
#1 Beam (5.5 cm)	8	#4 Beam (17 cm)	18
#2 Beam (8 cm)	8	#5 Beam (24 cm)	8
#3 Beam (11.5 cm)	18	Half Round Connector	14

When connected at 180° to a connector, two identical beams

two #1 beams connected at 180° have the same length as a #3

Half Round Screw Connector #5 Beam #4 Beam #3 Beam have a combined length equal to one longer beam. For example, #2 Beam #1 Beam #3 #3

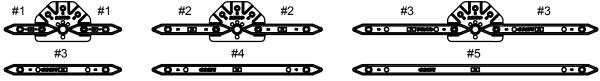

Truss Set Screws (ME-6994)

Each set of Truss Set Screws contains 75 Thumbscrews for attaching I-beams to connectors or load cells.

Cord Lock Spares (ME-6996)

A set of Cord Lock Spares includes 32 Cord Tensioning Clips (Cord Clips) and one roll of yellow braided cord. When attaching cords for lateral bracing or for suspension or cable-stayed bridges, Cord Clips are used to assist in adjusting the tension in the cords.

The Cord Clip does not come apart. It is best to thread the cord through the clip before the clip is installed on the bridge or structure. Prepare to thread the cord by holding the top half of the clip as shown in Figure A so the two



beam.

halves of the clip will separate, leaving an opening through which the cord is threaded. The cord is inserted into the end opposite the pointed end of the clip. The cord should be looped back through the clip as shown in Figure C. Then the Cord Clip can be used in the structure, using the attachment screw to tighten the clip shut. To adjust the cord tension, loosen the screw and pull on the cord to the desired tension and then tighten the screw.





Figure A: Hold half of the cord clip so the two halves separate

Figure B: Loop the cord back through the cord clip



Figure C: The cord goes around the screw hole



Figure D: The cord clip is ready to be attached to the structure using a screw

Full Round Connectors Spares (ME-6997)

Full Round Connector: The Full Round Connector has eleven slots, labeled A through H and X, Y, and Z, for attaching beams. There are six Full Round Connectors in the set.

Flat Round Connector: The Flat Connector has eight slots, labeled A through E, and X, Y, and Z, for attaching beams. There are six Flat Connectors in the set.


PAStrack Slot Connector: The PAStrack Slot Connector is a nut and bolt that allows a PAStrack to be connected to a structures model. There are six PAStrack Connectors in the set.


Full Round PAStrack Slot Connector


Axle Spares (ME-6998A)

The Axle Spares set includes two Axles each of three different lengths, twelve Pulleys, twelve O-rings, four Drive Wheels, four Tires, 24 Collets, and twelve Spacers.

The illustrations show example uses of axles, pulleys, and wheels.

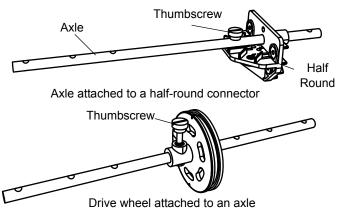

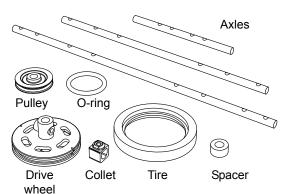


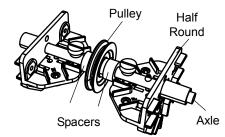


This pulley is free to rotate.

Angle Connectors (ME-6999A)

The Angle Connectors set includes 24 Straight Connectors, 24 Angle Connectors, and twelve Sliding Connectors. The Straight Connector can connect two beams to make a longer beam. The Angle Connector can allow a beam to be connected to a half round connector, full round connector, or flat round connector at an angle different than zero, 45, or 90 degrees. The angle connector also allows for a small adjustment of the length of the beam. The Sliding Connector allows one beam to be connected to another beam at any position along the length of the second beam..

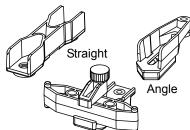

Sliding Connector on a #1 Beam

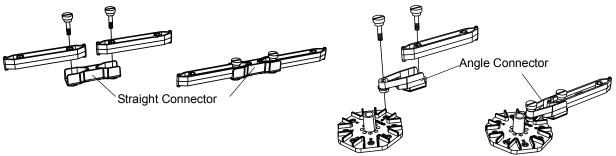

To use the sliding connector, loosen the thumbscrew and rotate the top "jaw" to the side. Place the beam onto the lower part of the connector, rotate the top "jaw" into place, and tighten the thumbscrew. The Structures Set includes two Angle Connector sets.

Connector Spares (ME-7002)

Each set of Connector Spares includes fourteen Half Round Connectors.

#6 I-Beam Spares (ME-7008)


The #6 I-Beam is 35 cm long. There are 24 beams per set.


Coaster Track (ME-9814)

The ME-9814 Coaster Track is a section of flexible track that is 9.1 m long. Side rails and center rails help to keep a Mini Car on the track.

Use a sharp knife or scissors to cut the track to the desired length.

Use a Track Coupler to connect two sections of Coaster Track together.

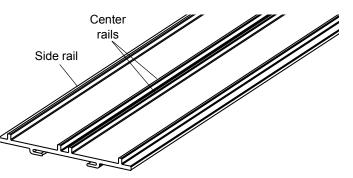

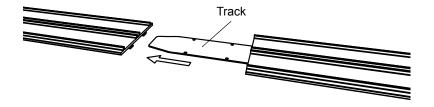

Slide the Track Coupler into the channel on the underside of each Coaster Track section.



Mini Cars (ME-9839 and ME-9840)

The ME-9839 Green Mini Car and ME-9840 Yellow Mini Car are designed to run on the Coaster Track.


If you are using a photogate (available separately) to measure the speed of the car, attach a photogate flag (included with the ME-6974) as pictured. The flag will interrupt the photogate twice as the car passes, allowing the speed to be calculated.


Mini Car Starter Bracket (ME-9856)


The ME-9856 Mini Car Starter Bracket is designed to help align the wheels of the Mini Car on the Coaster Track. Squeeze the metal clip of the starter bracket and insert it into the channel on the underside of the track.

Slide the bracket along the track to the desired position.

Drop the Mini Car into the bracket as shown.

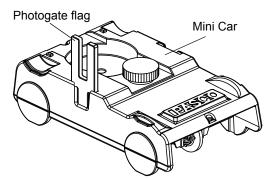

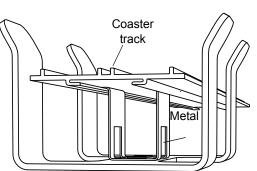

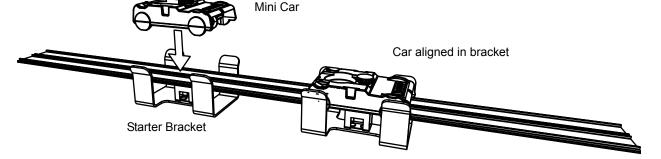

Total Quantities

Included Items	Qty	Included Items	Qty	Included Items	Qty
#6 Beam (35 cm long)	24	#4 Flexible Beam (17 cm)	18	Force Platform Structure	2
#5 Beam (24 cm long)	24	#3 Flexible Beam (11.5 cm)	18	Cord Tensioning Clip	32
#4 Beam (17 cm long)	54	Angle Connector	24	Yellow Cord	1 roll
#3 Beam (11.5 cm long)	54	Straight Connector	24	Coaster Track (9.1 m)	1
#2 Beam (8 cm long)	24	Sliding Connector	12	Mini Car (Green and Yellow)	2
#1 Beam (5.5 cm long)	24	Flat Round Connector	6	Mini Car Starter Bracket	1
Flat 2 X 3 Beam (12.5 cm)	16	Full Round Connector	6	Track Clip	24
Flat 3 X 4 Beam (19 cm)	16	Half Round Connector	70	Track Coupler	2
Flat #4 Beam (17 cm)	16	PAStrack Slot Connector	6	Mini Car Photogate Flat	2
#5 Flexible Beam (24 cm)	10	Screw (6-32)	450	Container (not shown)	2



Operation

Assembling Beams

All beams attach to connectors in the same way. Use the included screws (6-32, slotted) to attach beams to a connector (such as the half round connector) as illustrated.



Attaching beams to connectors


Adding Load Cells

To measure the compression and tension forces in individual members of a structure, add load cells (available separately) to the PASCO Structure. Replace a beam with two shorter beams and a load cell.

#5 beam = load cell + two #3 beams

#4 beam = load cell + two #2 beams

#3 beam = load cell + two #1 beams

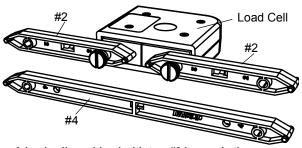

A load cell combined with two #2 beams is the same length as a #4 beam


Use thumbscrews to attach two beams to a load cell as shown in the figure.

When using load cells, assemble your structure with the screws loose. This will simplify the analysis by ensuring that the members experience only tension and compression without moments.

Example: Bridge with Load Cells

Static Load

The bridge shown in the figure incorporates six load cells to measure the tension or compression in various members. A hanging mass is used to apply load. The mass is adjusted so that the compression in one of the legs is 1.0 N. Compression is registered as a positive value and tension as a negative value.

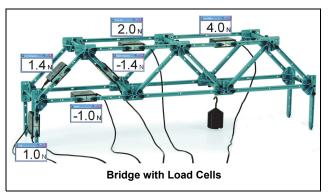

If the screws are loose, then the theoretical analysis

of the bridge can be carried out by assuming that the net force at each node is zero. Thus, the vertical component of compression in the left-most diagonal member must be 1 N (to oppose the force applied by the leg). The horizontal component must also be 1 N since the member is at a 45° angle. The predicted resultant force is:


$$\sqrt{(1.0 \text{ N})^2 + (1.0 \text{ N})^2} = 1.4 \text{ N}$$



Dynamic Load

With the load cells inserted as shown in the figure, push a Cart with its extra mass across the bridge. Zero the load cells before the measurement. Examine which members are under tension or compression.

Note that the Load Cell Amplifier can support up to six Load Cells in any combination (5 N, 100 N, or both).

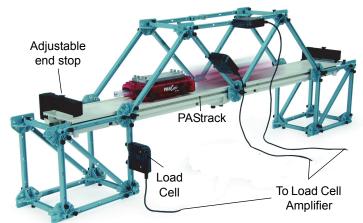


Recording the forces measured by the load cells as the cart traverses the PAStrack bridge.

Calibration.

See the instructions that came with the load cells for details about how to connect the load cells to an interface or datalogger. Follow the instructions to calibrate the load cells and use them to collect data. (Load cells are factory calibrated; however, you can re-calibrate them in software or on the datalogger. See the documentation for your software or datalogger for instructions.)

Large Structures Construction

The manual illustrates construction of twenty large structures such as trusses, bridges, and cranes. The descriptions show what components are used for each structure and how they are connected to each other. Close-up photographs show the details of construction where possible.

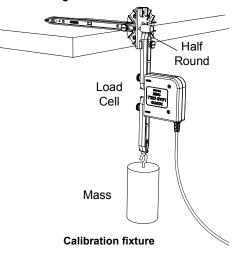


Structure	Structure	Structure	Structure
Car Building Jump	Suspension Bridge	Cantilever Truss	Arch Causeway
Cable Stayed Bridge	High Road/Low Road	Brachistochrone	Skyscraper
Arch Truss	Tied Arch Bridge	Drawbridge	House Frame
Roller Coaster	Baltimore Bridge	Vertical Lift Drawbridge	Angle Crane
Canyon Car Jump	Double Tied Arch Bridge	Tower Crane	Windmill

Tabl	^	-	- 1	
Iau	e			

Car Building Jump

Use beams, connectors, cord (for cross bracing), and a section of Coaster Track to build a ramp next to a building. Use Track Clips to attached the Coaster Track to the ramp. Add a Mini Car Starter Bracket to the top of the ramp and let a Mini Car roll down the ramp and "jump" onto the adjacent building's landing platform.

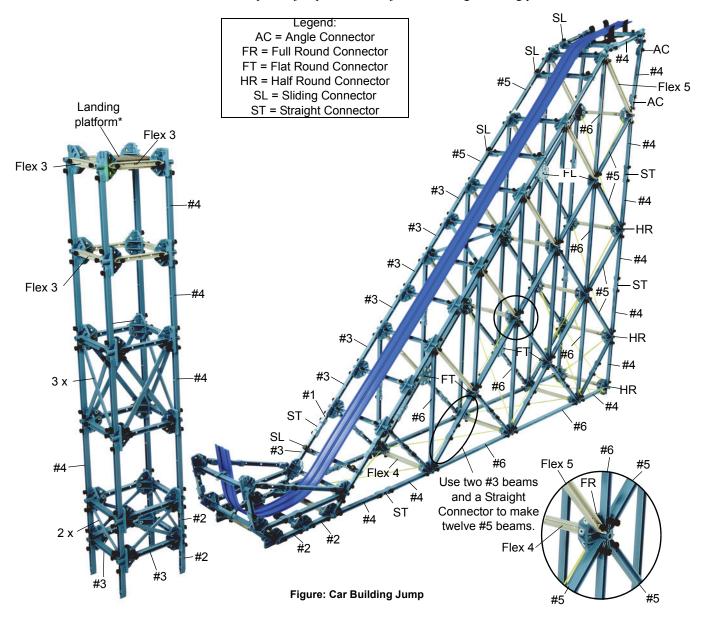


Adjust the position of the building as needed so that the car lands on top of the building.

*Make a landing platform using a square of cardboard or foam core board.

If the building is too high or too low, change the beams at the bottom of the building. Replace the #2 beams with #1 beams, or remove the bottom beams altogether.




Car Building Jump Details

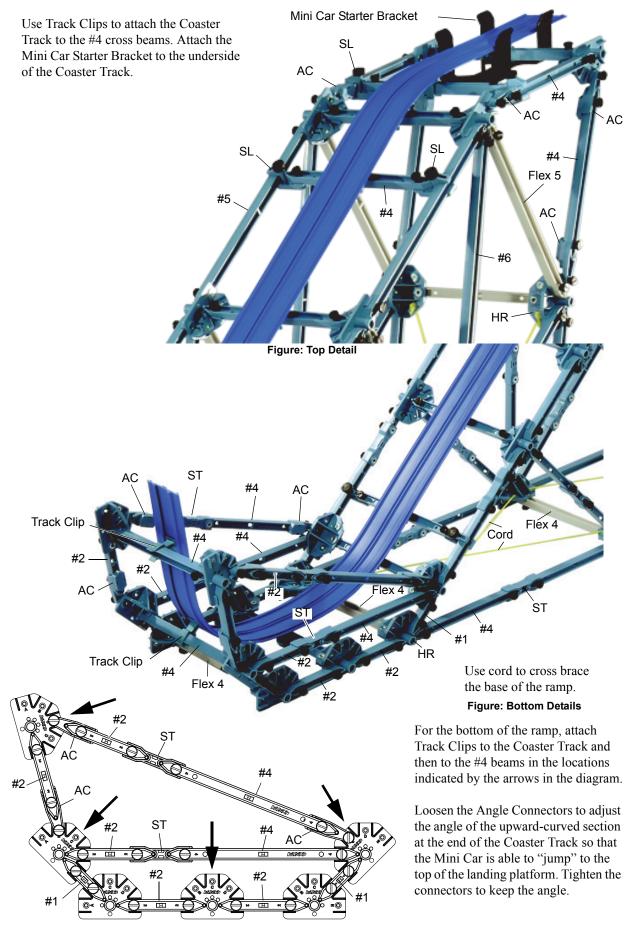

|245360®

Car Building Jump Details

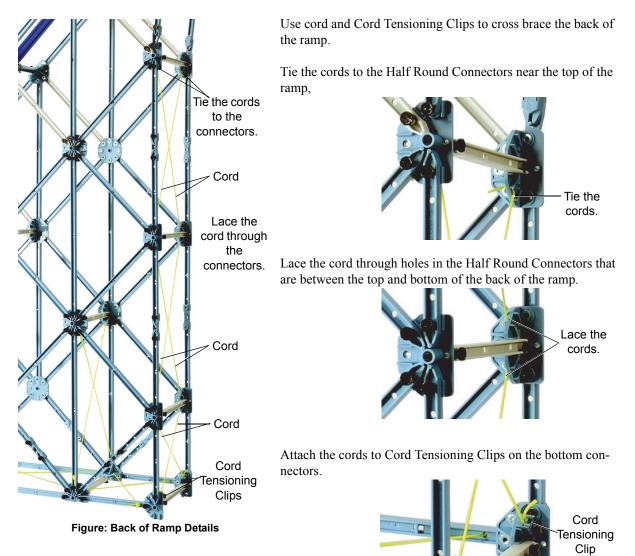

Cord Tensioning Clip



Cable Stayed Bridge

The Cable Stayed Bridge has a deck that is suspended on cords from two towers. There is a deck support structure at each end of the deck.

The modular approach to the construction of this structure is to build the deck in two halves and also build each tower and deck support structure before trying to suspend the deck. Do not attach the Coaster Track to the deck halves yet. Next, insert one half of the deck through one of the towers and use cord to suspend the deck from the top of the tower. Repeat for the other half of the deck. Connect the ends of the deck halves to the deck support structures and then connect the deck halves between the towers. Adjust the length of the cords so that the deck is level. Finally, use Track Clips to attach the Coaster Track to the finished deck..



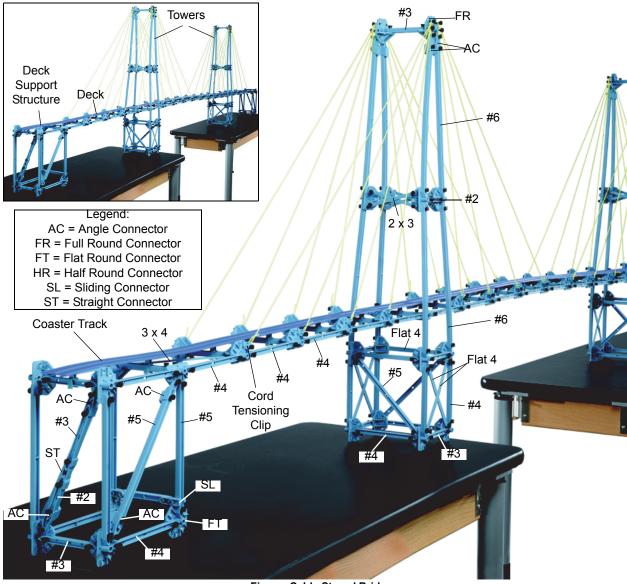


Figure: Cable Stayed Bridge


Cable Stayed Bridge Details

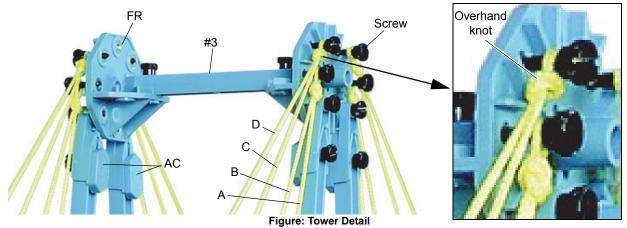

There are eight strands of cord from each side of the top of the tower that are used to suspend each half of the deck. Start with the strands of cord closest to the tower (for example, A and B in the figure). Get a piece of cord that is twice as long as the distance from the top of the tower to the deck. Make an overhand knot near the middle of the piece of cord and fasten the loop of the knot to the Full Round Connector at the top of the tower using a screw as shown. Put the two ends of the cord into Cord Tensioning Clips and attach the clips to two of the Half Round Connectors on the side of the deck. Repeat the process for strands C and D and the rest of the strands of cord.

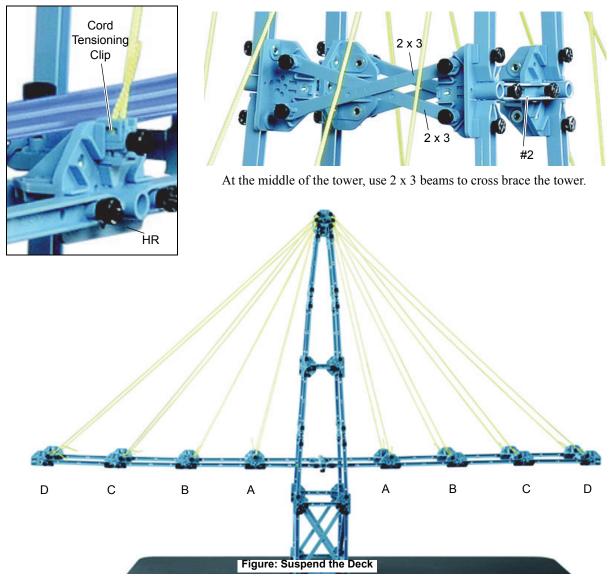


Arch Truss Bridge

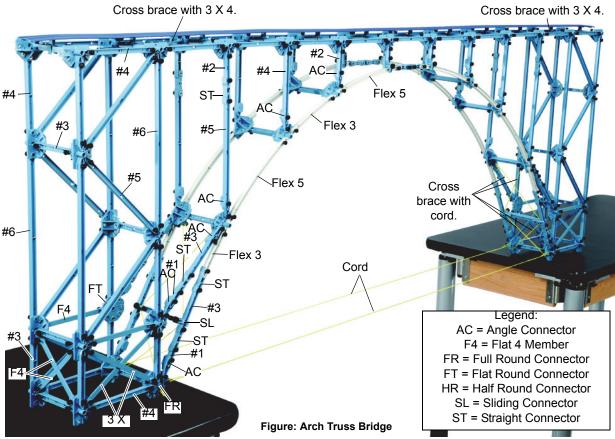

The Arch Truss design is one of the most frequently used designs in history. Construct the end supports and then add the deck. Tie cords from the base of one end support to the base of the other end support. Next, add the arch and suspension beams. Use Flat 3 X 4 members to cross brace the deck at the point where it joins the top of the end supports. Use cord to cross brace the bottom sections of the arch. Finally, use Track Clips to attach the Coaster Track.

Connect the center of the arch with Straight Connectors and a #1 Beam.


Use Sliding Connectors and #1 Beams to suspend the center of the arch.

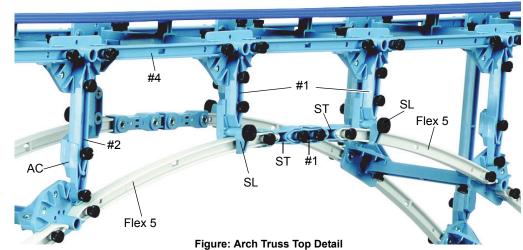


Arch Truss Bridge Details

Use two pieces of cord to cross brace the two lower sections of the arch. Attach the cords with Cord Tensioning Clips at one end. Lace the cords through the Sliding Connectors. Tie the cords to holes in the Full Round Connectors at the other end..

Repeat for the other side of the bridge.

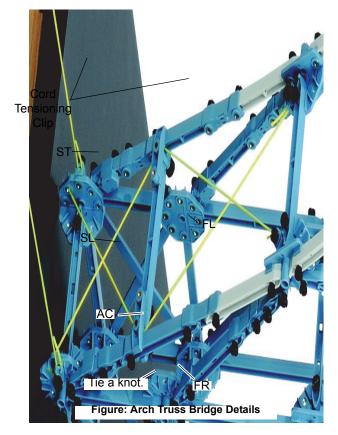




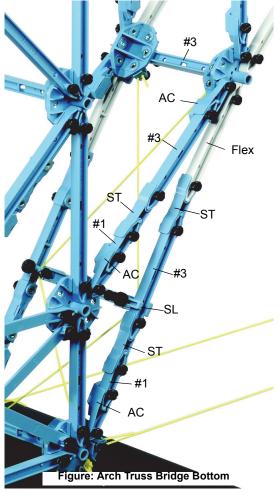


Roller Coaster

Use set components to build a "loop-the-loop" roller coaster model.

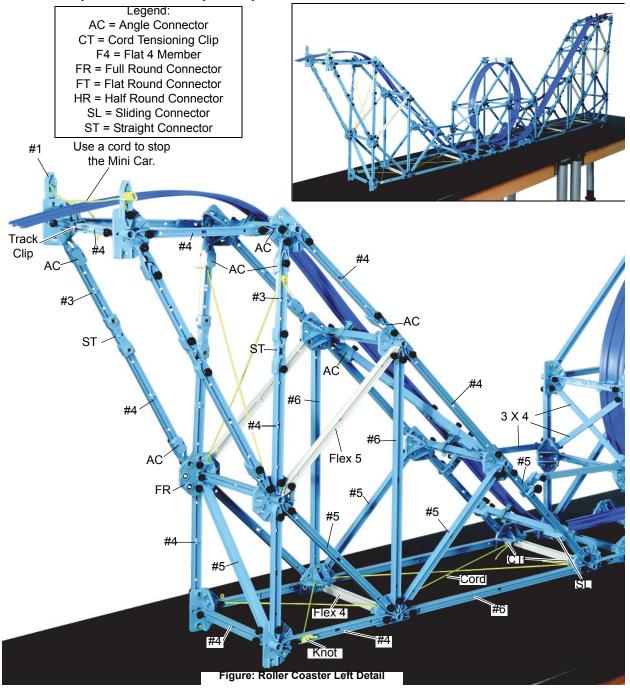

Roller Coaster Details - Right End

Make fourteen extra "#5" Beams using fourteen Straight Connectors and twenty-eight #3 Beams. Use cord to cross brace the bottom and the back of the ramp.

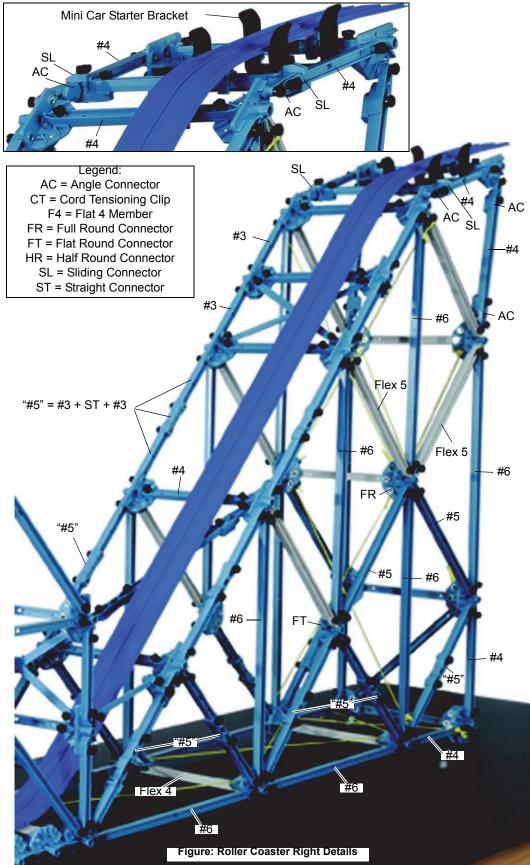


Roller Coaster Details - Center

Attach the Track Clip to a Flexible #4 Beam at the top of the Roller Coaster.



Figure: Roller Coaster Center Details

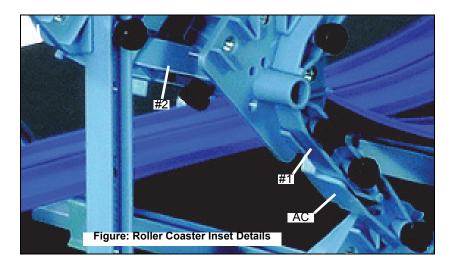

Canyon Car Jump

Use beams, connectors, clips and cord to construct a landing ramp.

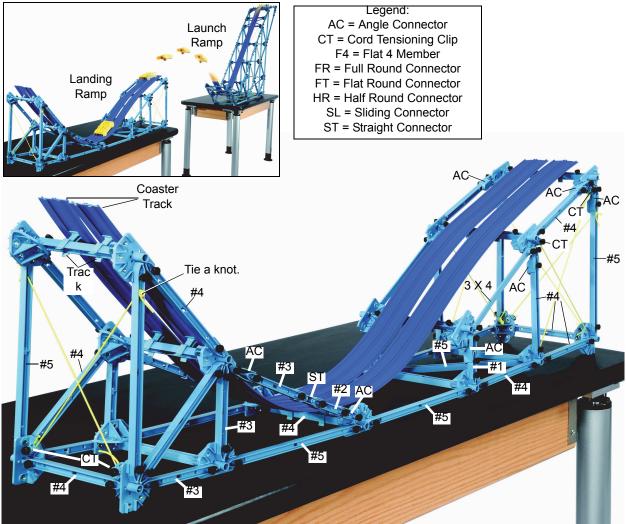


Figure: Canyon Car Jump Landing Ramp

Use cord and Cord Tensioning Clips to cross brace both ends of the landing ramp. Also use cord to cross brace the side panels at the right-hand end of the landing ramp. Use Track Clips to attach two sections of Coaster Track side-by-side to the #4 Beam cross members.

Canyon Car Jump Launch Ramp

Use connectors, cord, clips, and beams to construct a launch ramp. Include a Mini Car Starter Bracket at the top of the ramp

Make fourteen "#5" Beams using fourteen Straight Connectors and twenty-eight #3 Beams.

Use cord and Cord Tensioning Clips to cross brace the bottom and back of the launch ramp.

Use Track Clips to attach the Coaster Track to the #4 Beam cross members.

Add the Mini Car Starter Bracket.


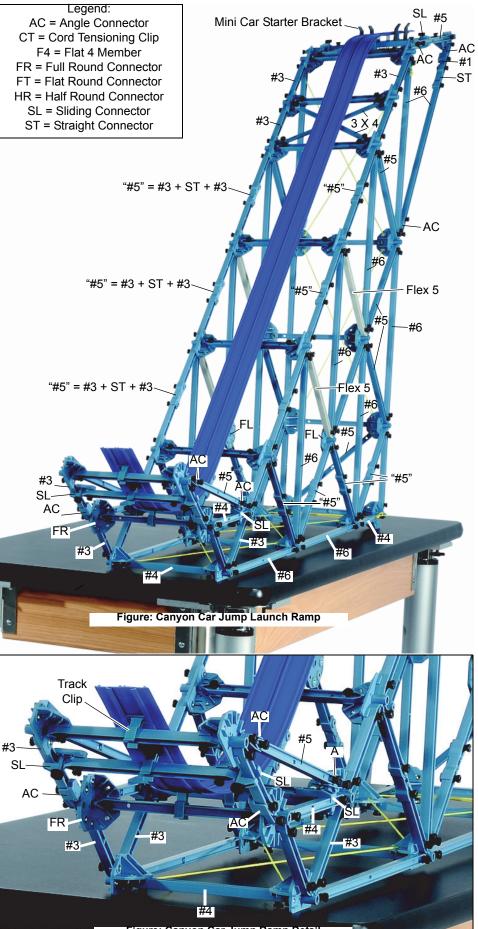


Figure: Canyon Car Jump Ramp Detail

Suspension Bridge

The construction of the Suspension Bridge uses a modular approach. The construction involves building the End Supports, Towers, Arches, and Deck, and then suspending the Deck from the Arches and connecting the Deck to the End Supports.



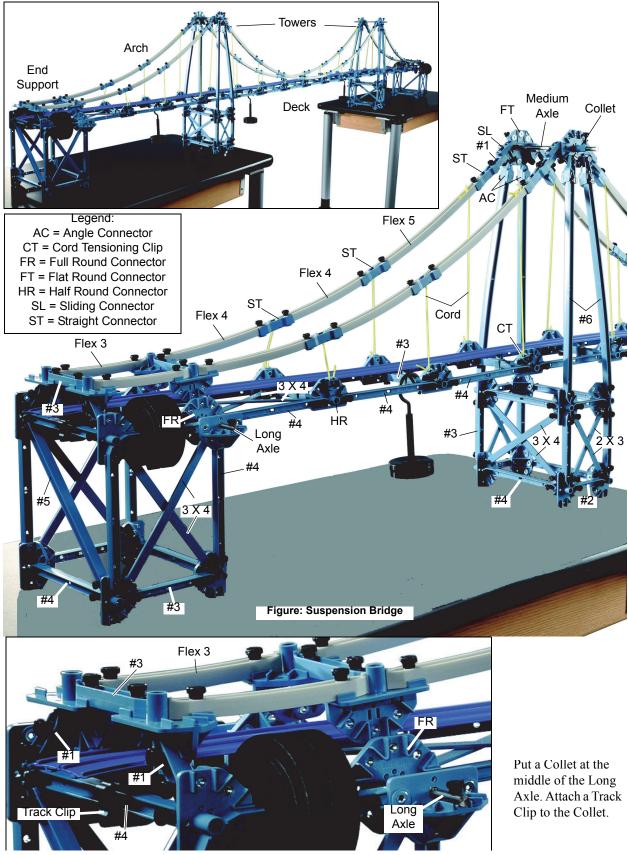

Suspension Bridge Details


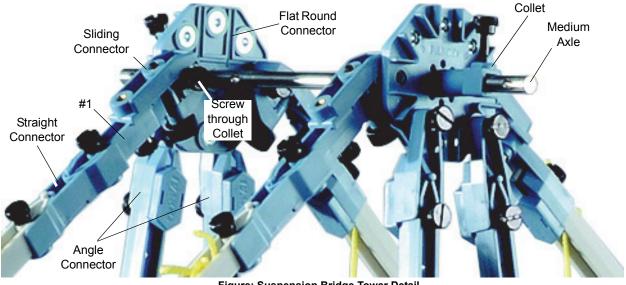



Figure: Suspension Bridge Tower Detail

Put a screw through the center hole of a Collet and attach the screw to the Flat Round Connector. Clamp the jaws of the Sliding Connector over the Collet. Connect a #1 Beam to the Sliding Connector and connect a Straight Connector to the #1 Beam. Repeat for the other arches.

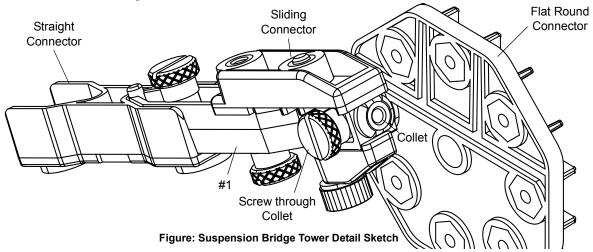


Suspend the deck from the arches using cord and Cord Tensioning Clips. Tie one end of each cord to the arch and attach the other end of the cord to the Cord Tensioning Clip.

Connect the end of the deck to the end support. Put one end of a Long Axle through the Half Round Connector of the end support and the Full Round Connector of the deck, Add a Collet to the Long Axle and then put the Long Axle through the Full Round Connector and Half Round Connector on the other side to attach the end of the deck to the end support. Position the Collet at the middle of the Long Axle and use a Track Clip to attach the Coaster Track to the Collet on the axle.

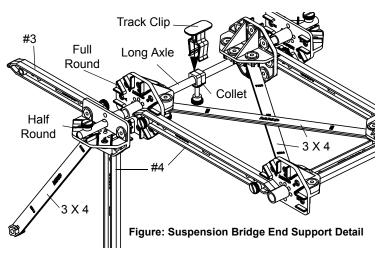


High Road/Low Road

Build a support for two Coaster Tracks side-by-side. One Coaster Track (the "High Road") has a constant gradual slope from start to finish. The other Coaster Track (the "Low Road") has a steep drop near the start and an equally steep rise near the end. If two Mini Cars start together on the two "roads", which car finishes first?

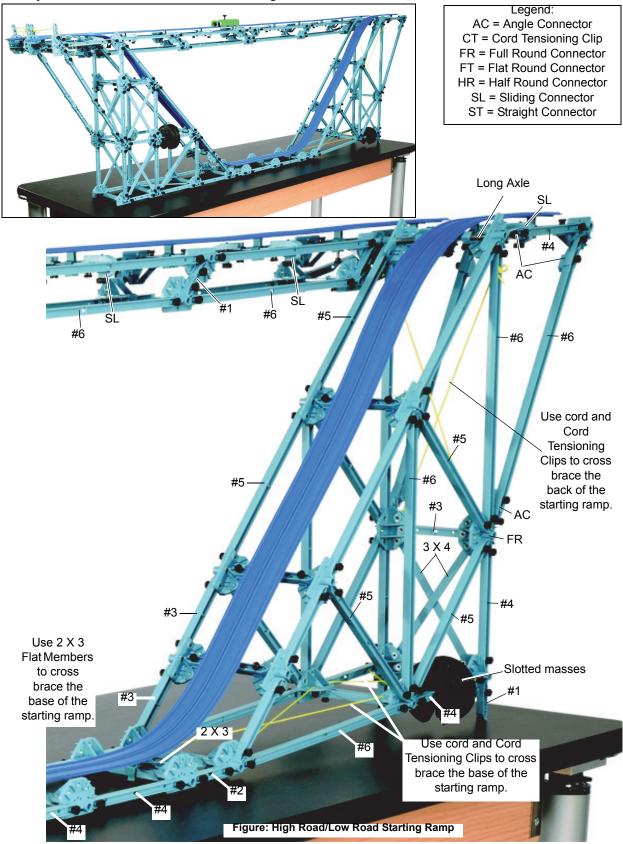


High Road/Low Road Details

The Long Axle extends through the support structures at the starting ramp for both the High Road and the Low Road.

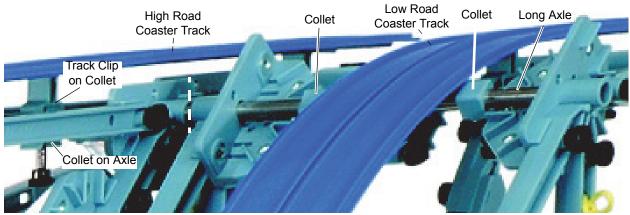

This illustration shows the High Road/Low Road starting ramp from another angle.

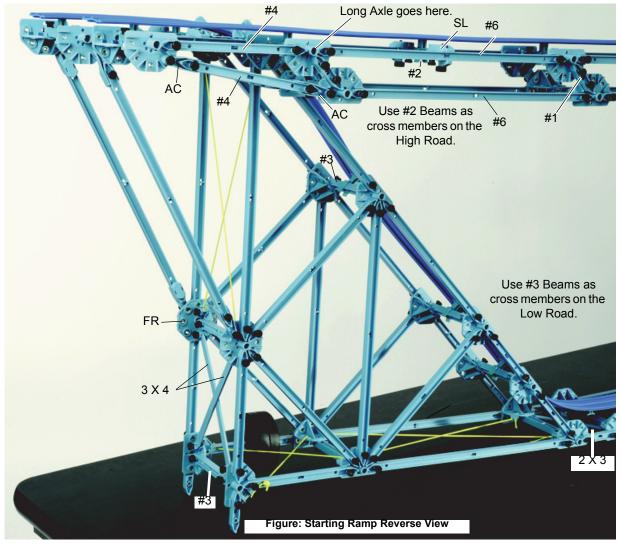

High Road/Low Road Details


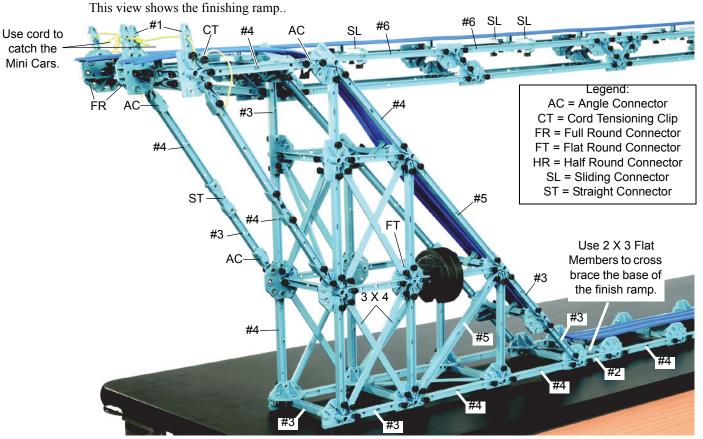


Figure: High Road/Low Road Finishing Ramp


Tied Arch Bridge



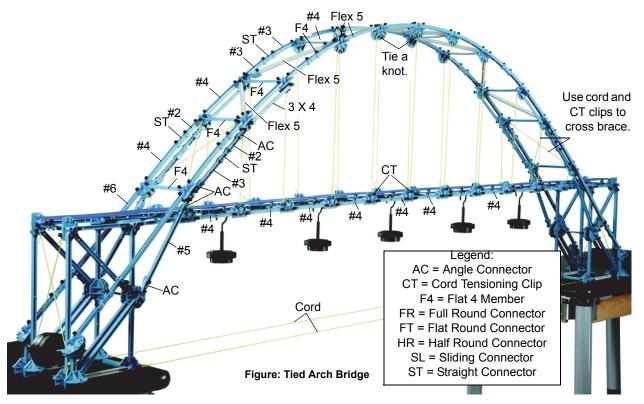


Tied Arch Bridge Details

Use a #3 Beam and two Sliding Connectors to support the slotted masses at the back of the end support.

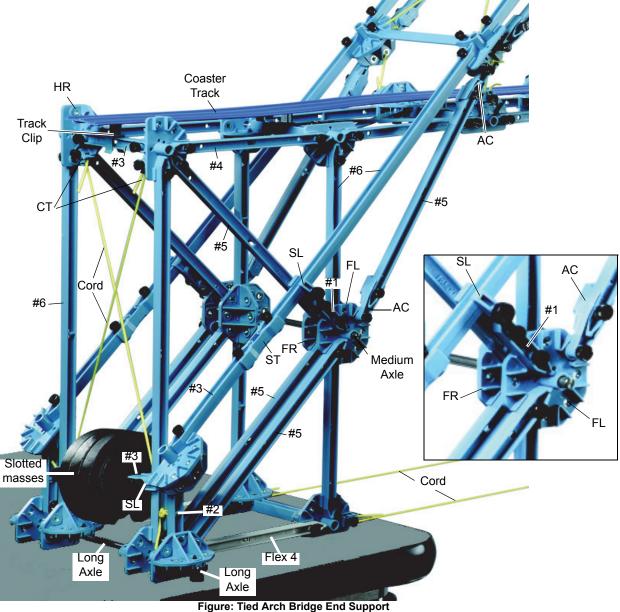


Use cord and Cord Tensioning Clips to cross brace the back of the end support. Tie cords between the bottoms of the two end supports.

Tied Arch Bridge Details

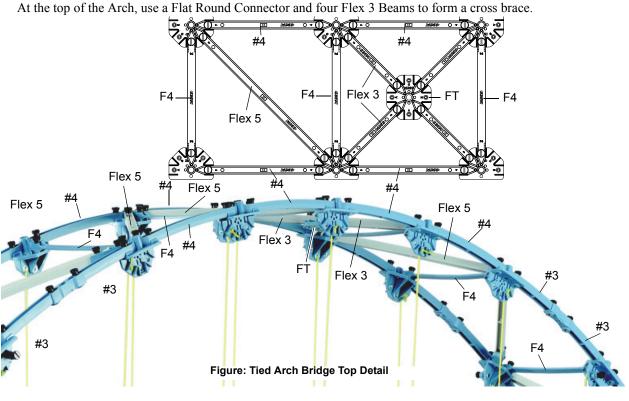

Baltimore Bridge

Use 40 #3 Beams and twenty Straight Connectors to make twenty extra "#5 Beams".

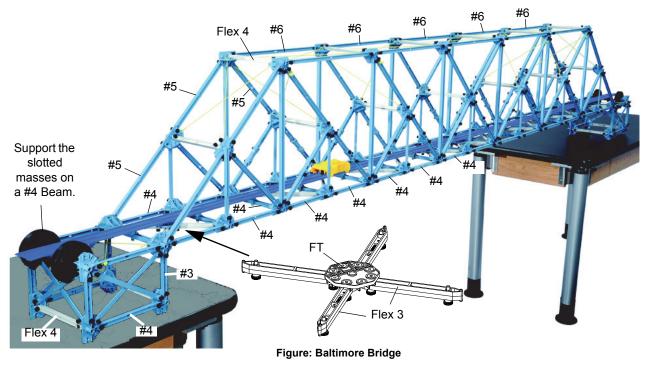


Use Flat Round Connectors and four Flexible #3 Beams to make four cross braces for the deck. Place one at each end of the deck where it joins the end support and put the other two on both sides of the center of the deck.



Baltimore Bridge Details

Use cord and Cord Tensioning Clips to add cross bracing to each panel on the top of the truss and to several of the diagonal panels. Use Flexible #5 Beams to cross brace several of the panels of the deck.

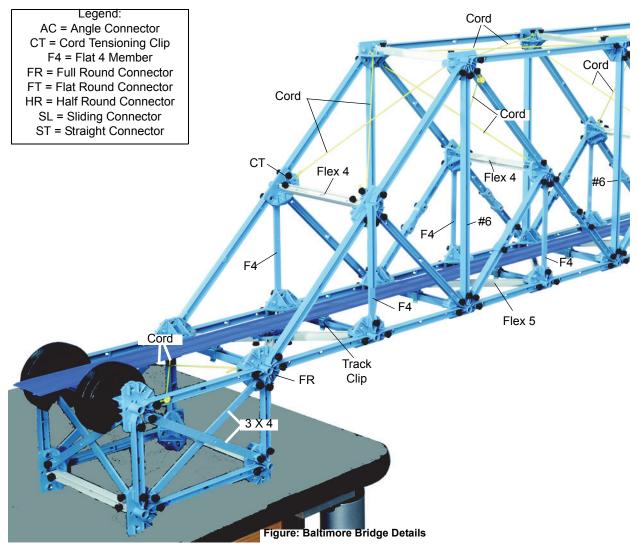

Baltimore Bridge Details

Use a Full Round Connector at the center of the bridge.

Add two of the cross brace structures to the deck on both sides of the center of the deck.


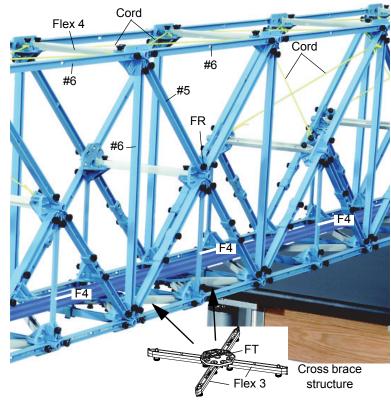


Figure: Baltimore Bridge Details



Double Tied Arch Bridge

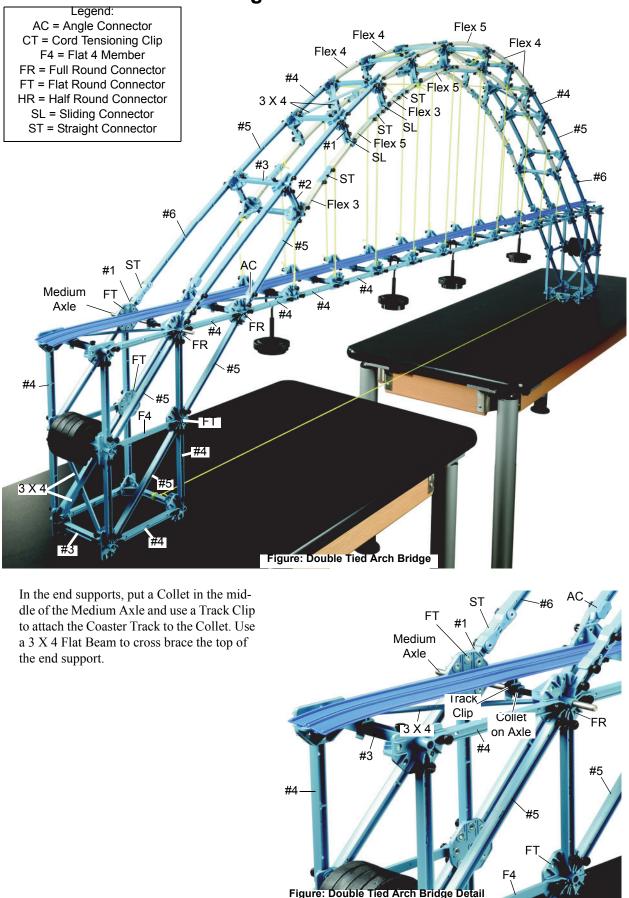

Double Tied Arch Bridge Details

Flex 5 Flex 4 Flex 4 Flex 4 Flex 4 3 X 4 Flex 5 #/ SL #4 ST Flex 4 Flex 3 SL ST Flex 5 Figure: Arch Details SL Note that there are three different positions on the Half Round Connector for the Cord Clip. Choose Cord the position that makes the cord as vertical as possible.Adjust the length of the cords to make the deck level. Add Masses СТ The mass hanger usually hangs from the #3 Beam CT that is near the #1 Beam where the 5 N Load Cell will be attached.Re-adjust the tension in the supporting cords to make the Deck level and give the bridge is proper shape.


Use cord and Cord Clips to support the deck from both sides of the arch..





Cantilever Truss Bridge


The Cantilever Truss Bridge balances the torque on one side of the center tower with the torque on the other side.

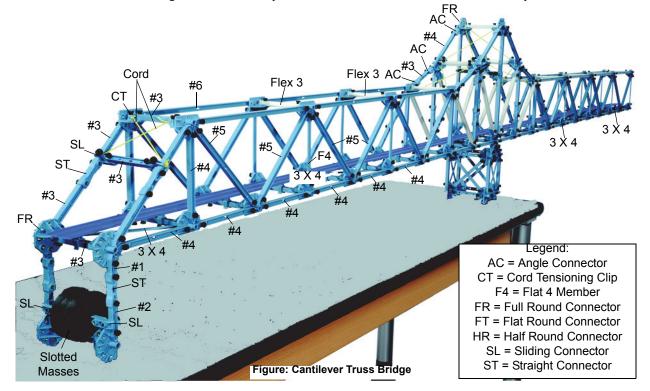

Use cord and Cross Tensioning Clips to cross brace the top section of the truss at the end of the bridge. Use Track Clips to attach the Coaster Track to the #3 Beam cross members. On each "arm" of the bridge, use a Flat 3 X 4 Member to cross brace the first section and the fifth section (counting from the end toward the tower).

Cantilever Truss Bridge Tower Details

Use cord and Cross Tensioning Clips to cross brace the top sections of the truss tower. Use Full Round Connectors at the top and the bottom of the #6 Beams at the center of the tower. Use two Sliding Connectors and a #3 Beam as a cross member between the #6 Beams at the center of the tower.

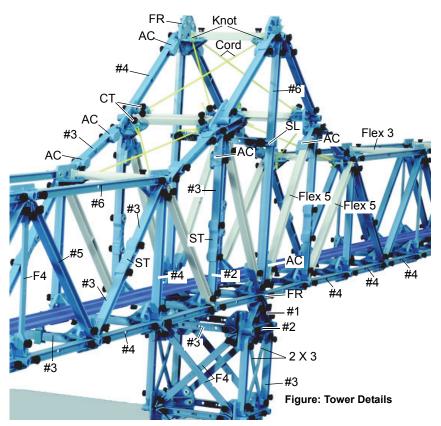


***/_{_**®

Brachistochrone

The word "brachistochrone" is from the Greek words meaning "shortest" ($B\rho\alpha\chi\iota\sigma\tau\sigma\zeta$) and "time" ($\chi\rho\sigma\nu\sigma\zeta$).

Imagine two fixed points in a vertical plane so that one of the points is not directly below the other. If a particle starts from rest at one of the points and descends to the other point under its own weight, the *brachistochrone* is the path the particle must take to reach its destination in the briefest time. The *brachistochrone* may not be the shortest path, but it is the quickest path.

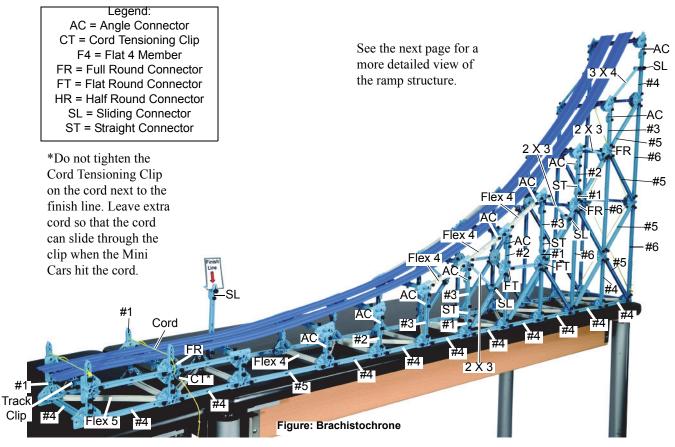


Set up the structure with two Coaster Tracks side-by-side. Place two Mini Cars at different starting points on the track and let them go. What happens?

Use cord and Cord Tensioning Clips to cross brace the back of the high end of the ramp.


The brachistochrone curve is part of an inverted cycloid. A cycloid is the path of a point that is fixed on the edge of a circle as the circle rolls along a straight line.

Imagine a straight ramp and a brachistochrone curve as shown. In a race from A to B between a ball rolling down the brachistochrone curve and a ball rolling down the straight ramp, which ball will finish first?

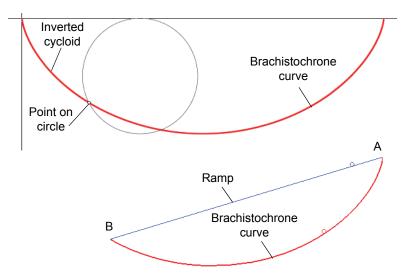




Brachistochrone Details

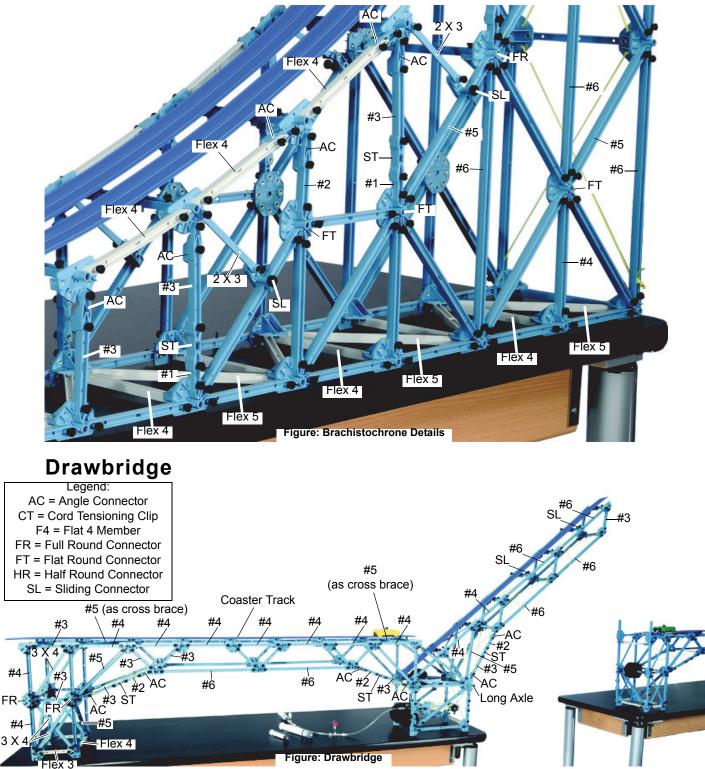


The drawbridge structure is a "working" model. You can use parts from the ME-6984 Structures Hydraulic System (not included*) in order to raise and lower the drawbridge. Use #4 Beams or Flexible #4 Beams as cross members on the bridge support section. Use two #3 Beams and a Straight Connector to build extra #5 Beams. Use two #2 Beams and a Straight Connector to make extra #4 Beams. Note: Additional members from PASCO Structures Sets are needed in order to extend the right hand end of the bridge shown in the illustration.

(*See the PASCO catalog or the PASCO Web site at www.pasco.com for information about the ME-6984 Structures Hydraulics System.)

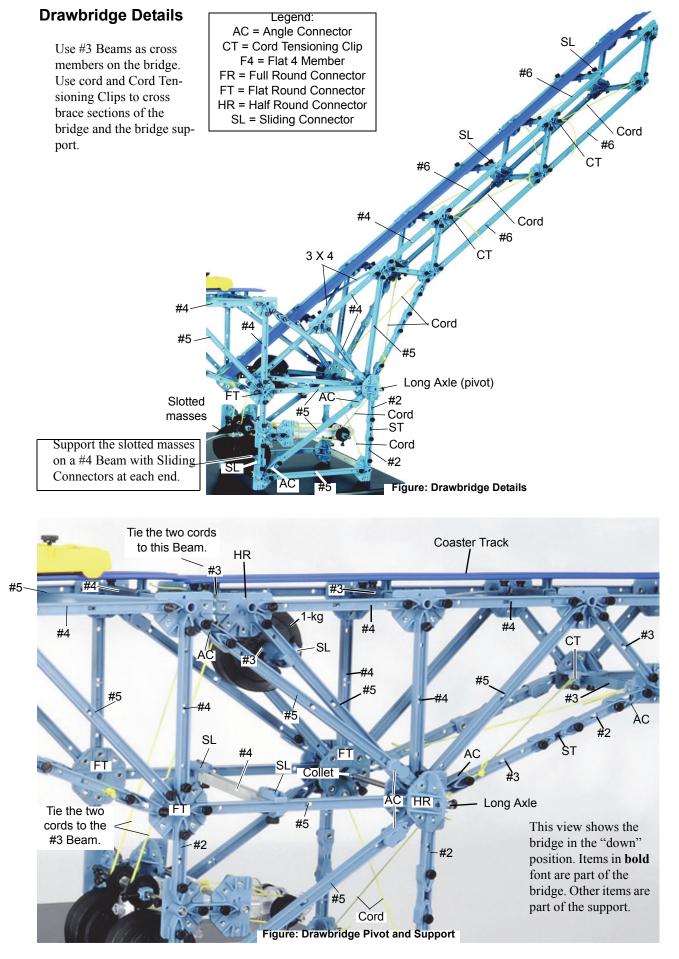



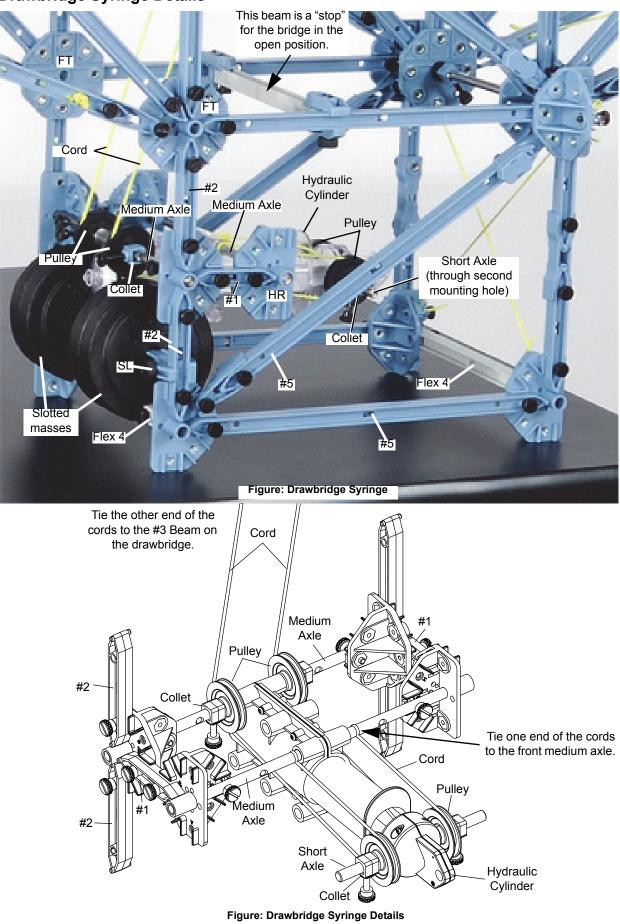

Drawbridge Details Continued

Use a #4 Beam or Flexible #4 Beam and two Sliding Connectors to form a "stop" bar for the drawbridge when it is in the closed position.



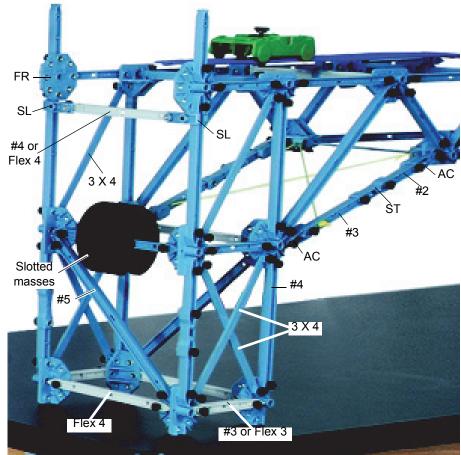


Figure: Drawbridge Right Side


Vertical Lift Drawbridge

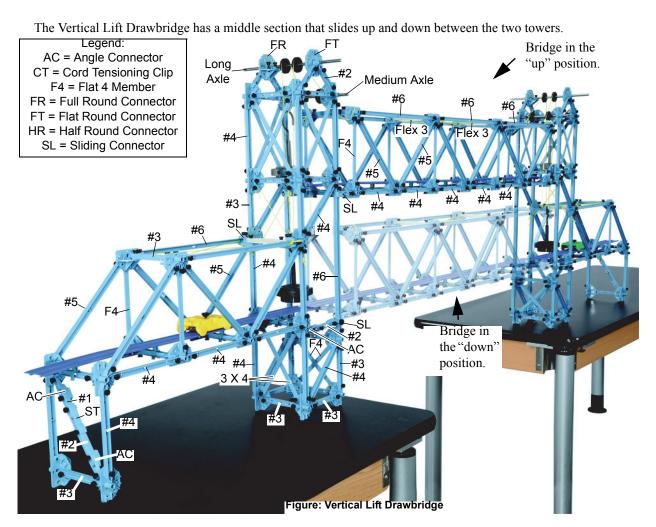

The middle section is counterbalanced by hanging masses connected by cords over pulleys. The hanging masses move up and down inside each tower.

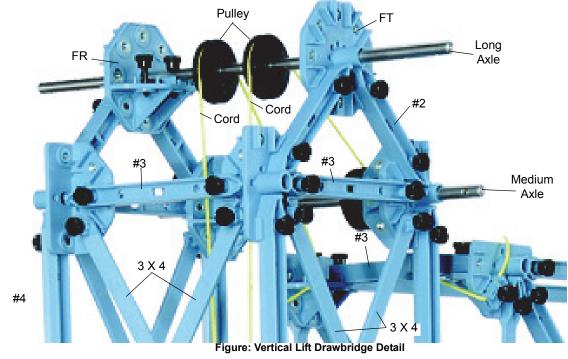

Vertical Lift Drawbridge Details

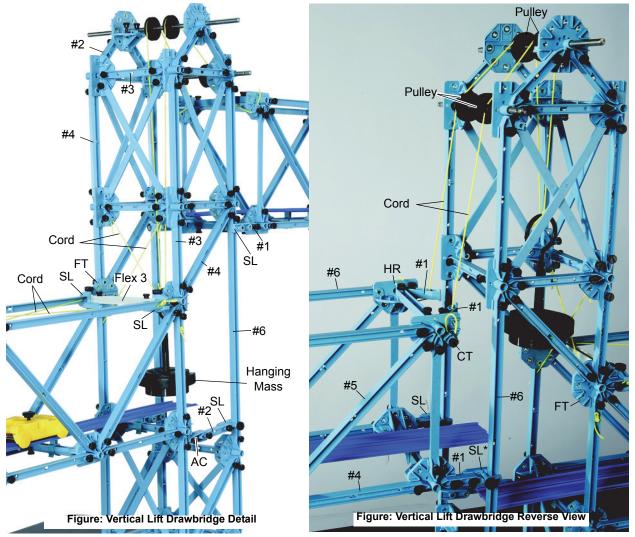


(*Do not overtighten the Sliding Connectors that grasp the vertical #6 Beams. The middle section must be able to slide up and down the #6 Beam.)

Use cord and Cord Tensioning Clips to cross brace the tower and the tower support. For the hanging mass cord, determine how long the cord must be so that the middle section can slide up and down on the #6 Beam on the outside of the tower while the hanging mass moves up and down inside the tower. Double that length. Thread the ends of the cord through the #1 Beams at the top edge of the middle section. Fasten the ends of the cords to Cord Tensioning Clips on the Half Round Connectors. Place the cords over the pulley, and hang the mass hanger from the middle of the cord. Adjust the amount of mass on the mass hangers so that the middle section of the bridge can be lifted or lowered easily.




Tower Crane


The Tower Crane stands 2.3 meters tall and can be mounted on a PASPORT Force Platform (PS-2141 or PS-2142) or a ScienceWorkshop Force Platform (CI-6461) using the Force Platform Structure Brackets (ME-6988).

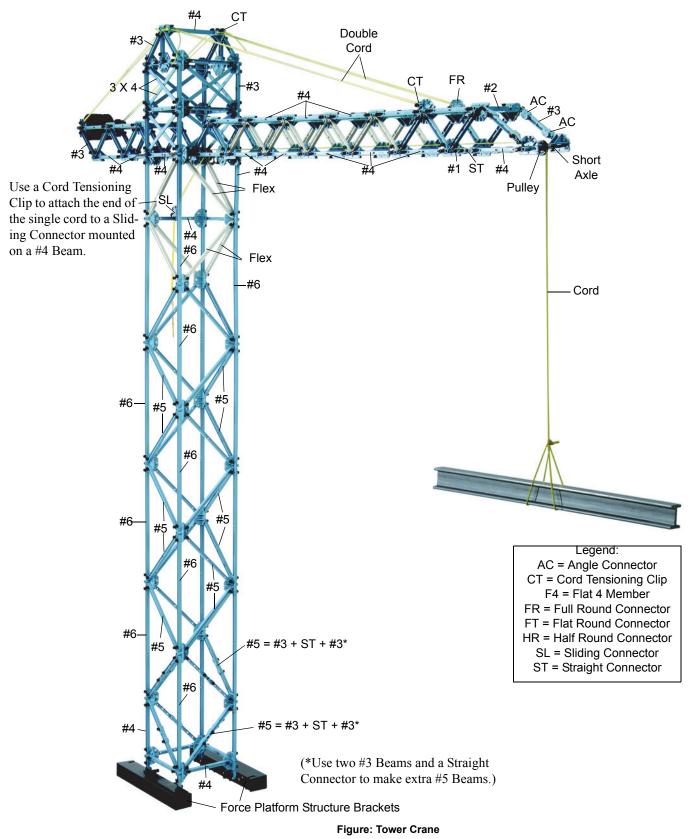

The Tower Crane consists of three parts: tower, rear boom and front boom. The tower has a #4 Beam by #4 Beam cross section, uses #4 and #6 Beams for the vertical members, and #5 Beams for the cross braces. The rear boom is a #3 Beam in width and the front boom is a #2 Beam in width. Flexible beams are used when there are no more #5 or #3 Beams.

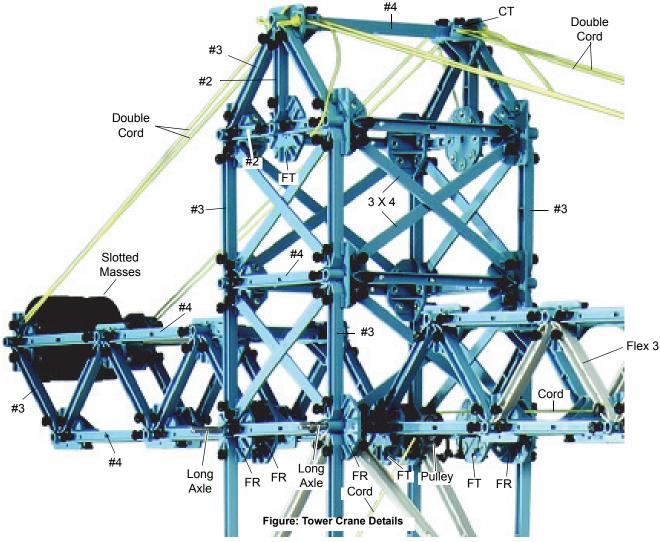


Put a Collet on each side of the Pulley to keep it in place on the front Long Axle. Make sure that the front Full Round Connectors are oriented so that the "fin" is vertical.

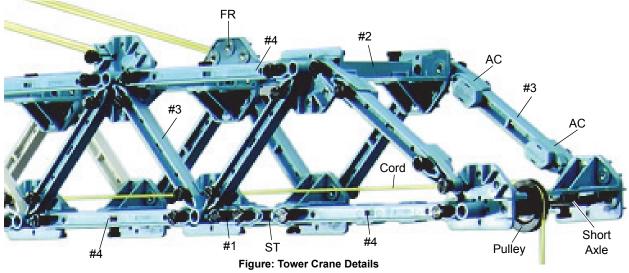




Tower Crane Boom Details

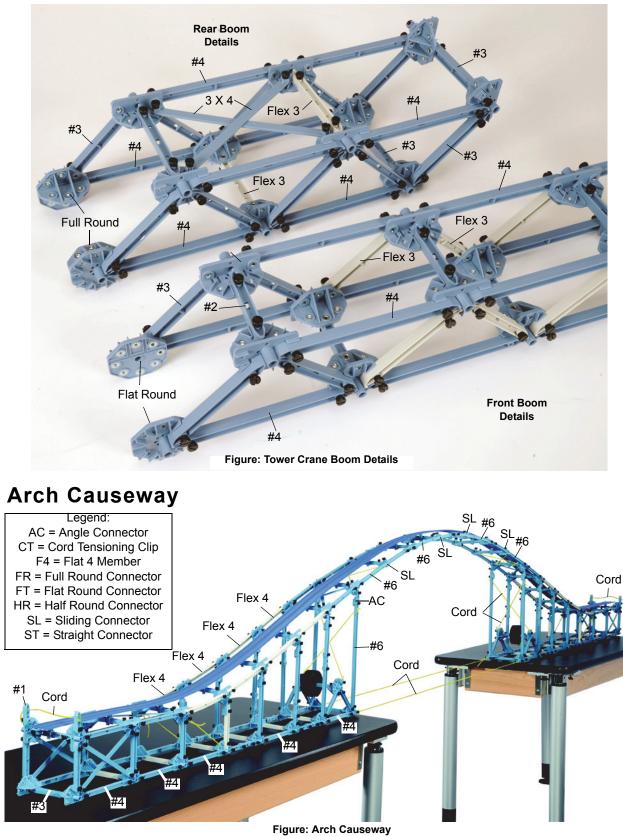

Use cord to connect the two support structures and to cross brace the tallest vertical sections. The arch uses four #6 Beams on each side with #3 Beams and Sliding Connectors as cross members.

Arch Causeway Details

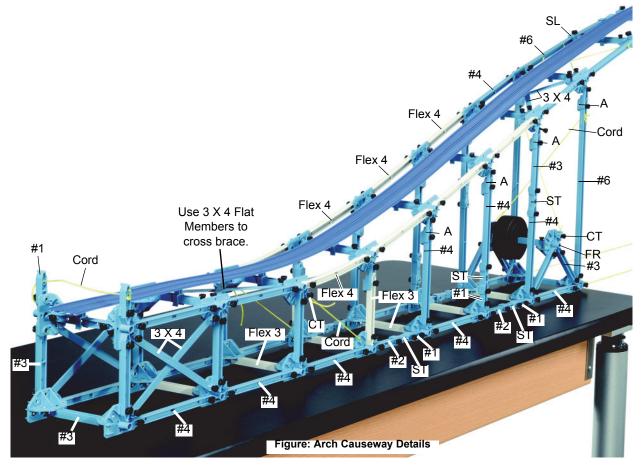

Use Flexible #3 Beams as cross members in the base of the support when there are no more #3 Beams.




Skyscraper

After the invention of the elevator, "skyscraper" buildings became a reality. Use #2 Beams to attach the bottom of the structure to the Force Platform Structures Bracket as shown.


The skyscraper structure models a sixteen story building. Use #4 Beams as the vertical members for the bottom eight stories and use #3 Beams as the vertical members for the top eight stories. Use #6 Beams and #6 Beams for the horizontal members of each story. Use two #3 Beams and a Straight Connector to make extra #5 Beams. Use two #4 Beams and a Straight Connector to make extra #6 Beams.

The skyscraper structure can be used to demonstrate what happens to a building during the shaking caused by an earthquake.

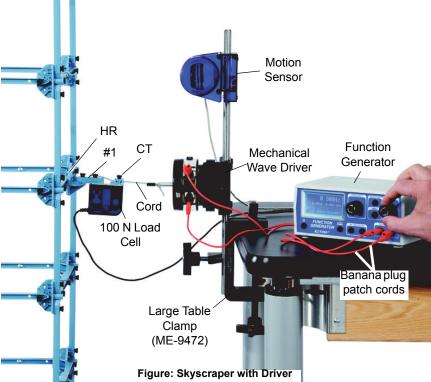


For example, replace a Straight Connector with a Half Round Connector. Add a #1 Beam to the Half Round. Attach a 100 N Load Cell (PS-2200) to the #1 Beam. Use cord to connect the Load Cell to the shaft of a Mechanical Wave Driver (SF-9324). Connect banana plug patch cords (SE-9750) from the wave driver to a Function Generator (PI-8127) and use the function generator to power and control the wave driver.

Mount a Motion Sensor (PS-2103) on a support rod (ME-8988). Connect the Motion Sensor and the 100 N Load Cell to a PASCO interface to record, display, and analyze the data as the skyscraper shakes.

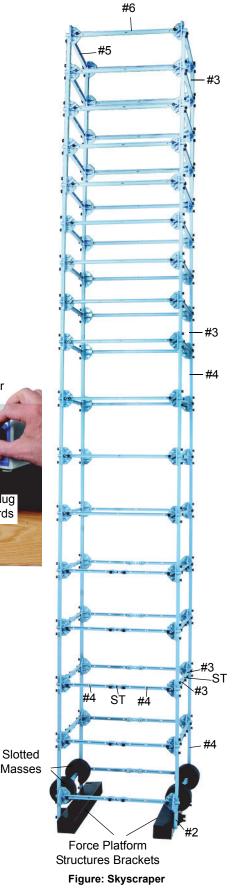


House Frame

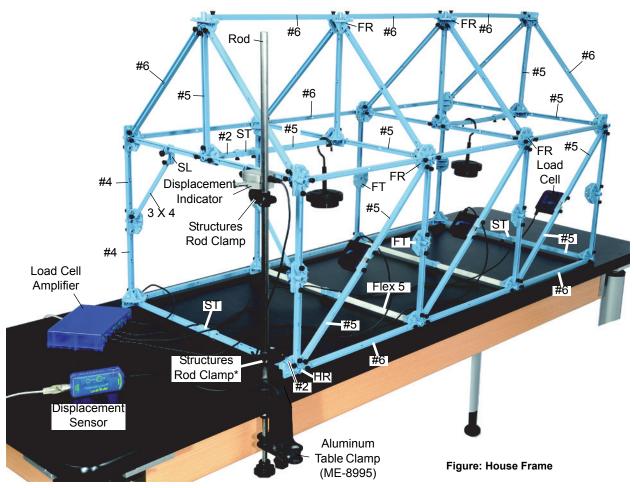

To hold the House Frame in place, attach a #2 Beam to the Half Round Connector at one corner. Put a clamp on the edge of the table and use it to hold a support rod (ME-8995). Attach a Structures Rod Clamp between the support rod and the #2 Beam at the corner of the House Frame.


To make measurements of the flexing of the House Frame when masses are hung from the members, mount the Displacement Indicator of the Displacement Sensor (PS-2204) on another Structures Rod Clamp and mount the clamp on the support rod so that the plunger of the indicator contacts a beam on the House Frame. To make measurements of the forces in the members, mount 100 N Load Cells as shown and connect them to a Load Cell Amplifier (PS-2198).

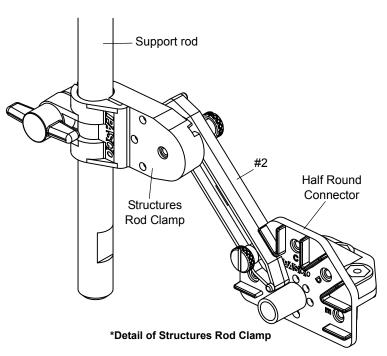





Angle Crane

The Angle Crane is one of the largest structures to be built with the Large Structures Set. The crane's boom can be raised or lowered using the Hydraulic Cylinder and other parts from the ME-6984 Structures Hydraulics System (available separately). The Angle Crane has three main parts: Boom, Tower, and Base. The Tower is joined to the Base by an axle mounted on the Base, and the Boom pivots on another axle near the top of the Tower.

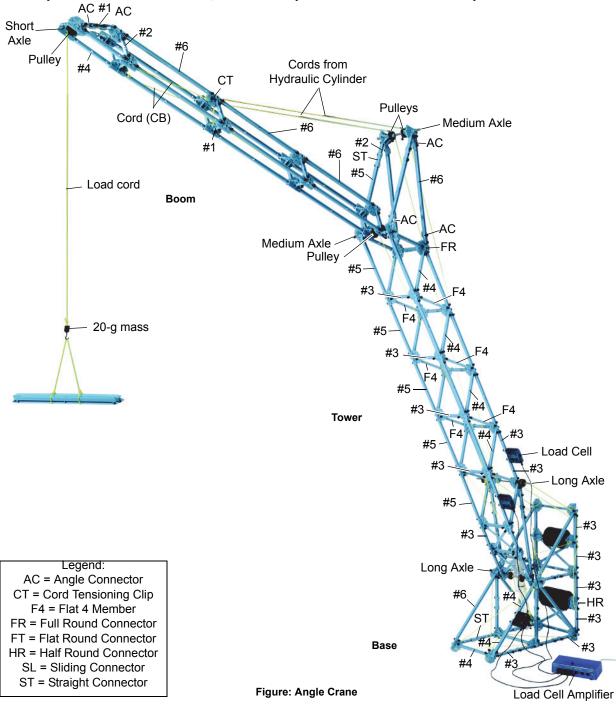


Use two #3 Beams and a Straight Connector to make ten extra #5 Beams for the Base and Tower. The hook on the end of the cord is a 20 g mass from the Hooked Mass Set (SE-8759). Slotted masses on the back of the base are counterweights for the Tower and Boom, and forces in the Tower and Base are measured by 100 N Load Cells connected to a Load Cell Amplifier. The Base width is a #4 Beam, the Tower width is a #3 Beam, and the Boom width is a #1 Beam.

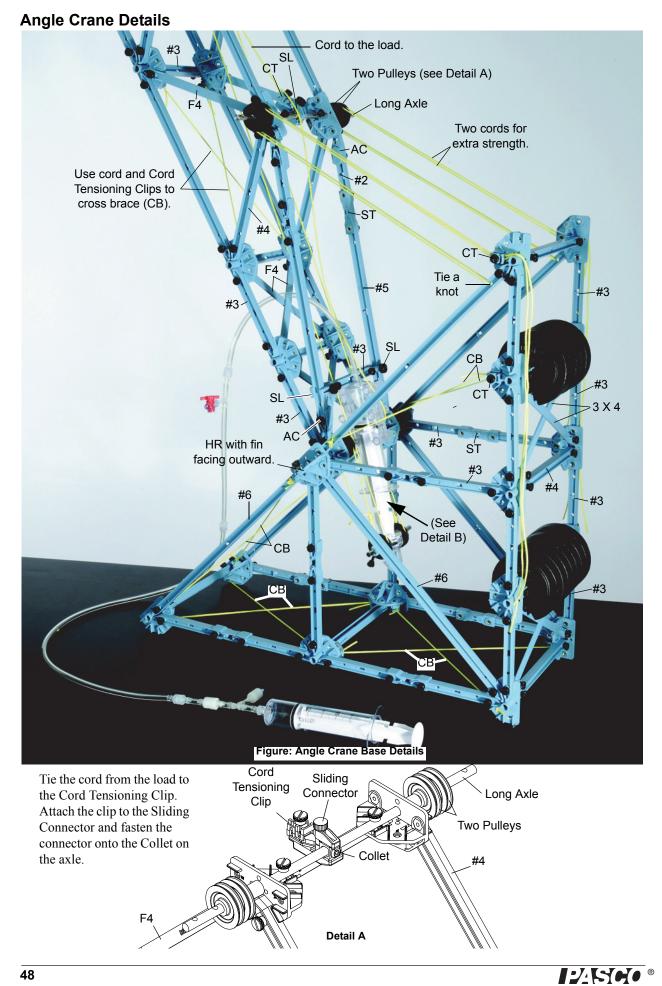

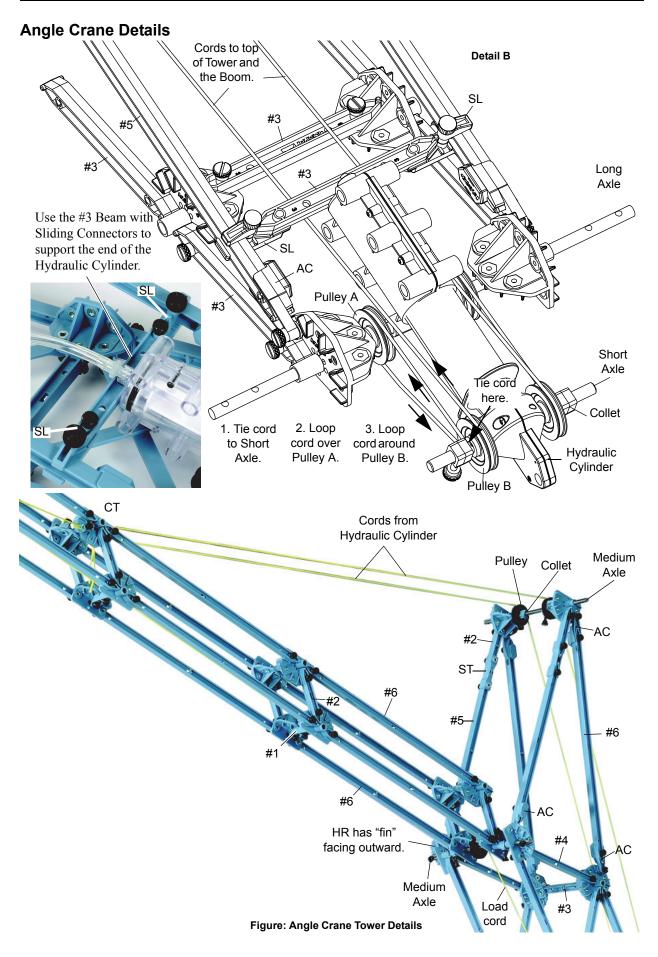

Windmill


The Windmill is another working model that can demonstrate how the kinetic energy of the wind can be converted into other forms of energy, including gravitational potential energy. In this case, the windmill turns an axle that winds up a cord attached to a hanging mass.

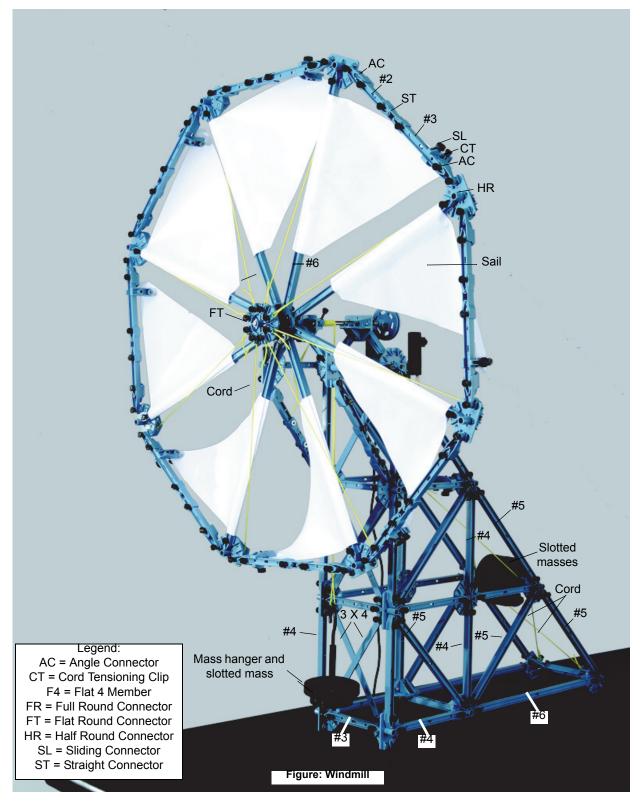

A template for the paper sails is included at the end of the instructions. Copy (or trace) the template and cut out the sail from a sheet of 8.5 by 11 inch (21.5 by 28 cm) paper.




Windmill Details

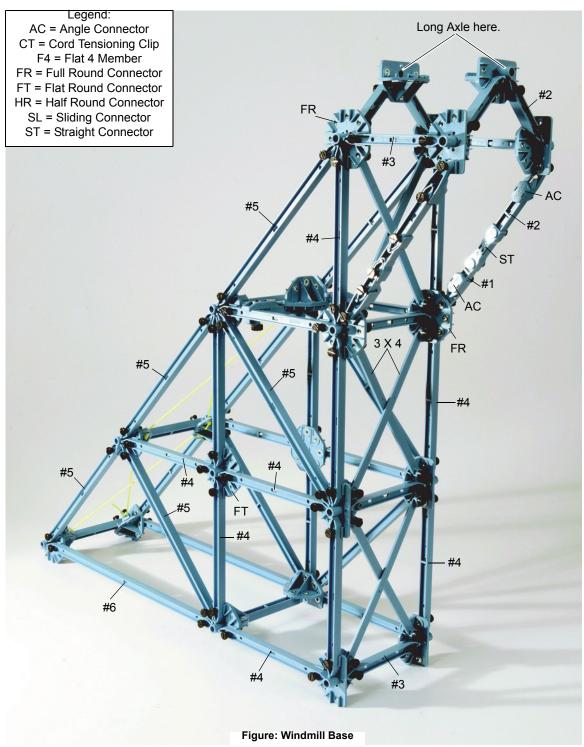


Make a sleeve in each sail to slide over the #6 Beams by folding the lower part of the sail to the dotted line, and then taping the edge in place. Clamp the small end of the sail between a Sliding Connector and a Cord Tensioning Clip.

Windmill Details

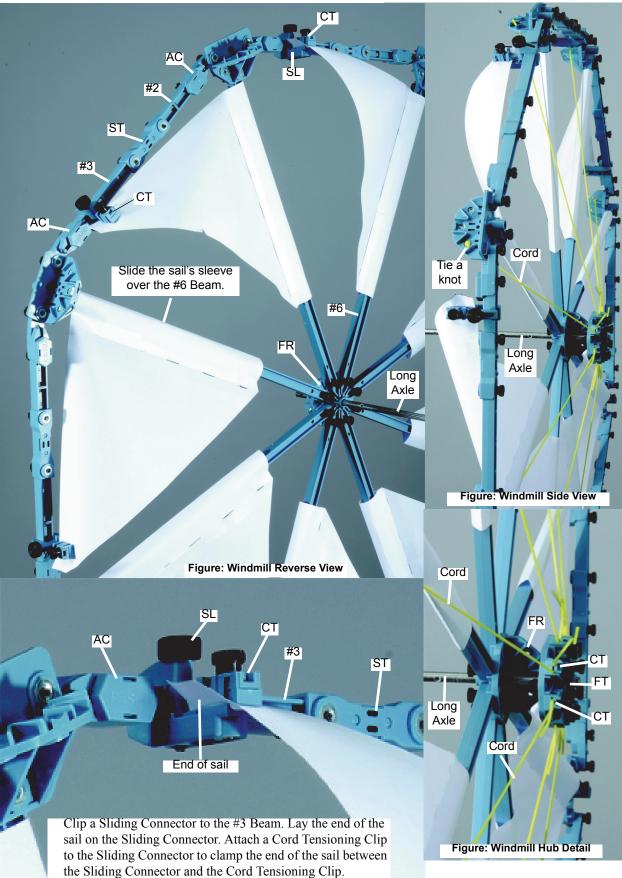

Windmill Details

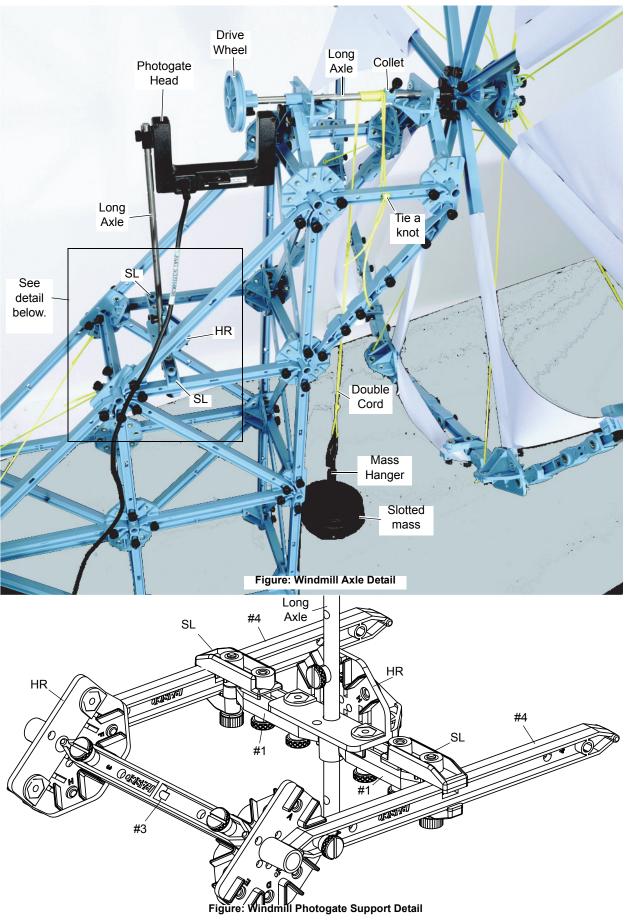

Technical Support

Before you call PASCO technical support, have the apparatus and this user's guide available. Please note the following:

- Product name and model number (e.g., Large Structures Set, ME-7003)
- Approximate age of the product;
- Detailed description of the problem/sequence of events required to duplicate the problem.

For assistance with any PASCO product, contact PASCO at:

Address:	PASCO scientific 10101 Foothills Blvd. Roseville, CA 95747-7100
Phone:	916-462-8384 (worldwide) 877-373-0300 (U.S.)
Web:	www.pasco.com
Email:	support@pasco.com

For more information about the Large Structures Set and the latest revision of this Instruction Manual, visit the PASCO web site and enter ME-7003 into the Search window.

Limited Warranty For a description of the product warranty, see the PASCO catalog.

Copyright The PASCO scientific 012-12670A *Large Structures Set Instruction Manual* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Trademarks PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.

Patents Pending: The following are some of the PASCO products that have patents pending:

ME-6990 Truss Set	ME-6991 Bridge Set
ME-6992A Advanced Structures Set	ME-6993 Truss Set Members
ME-6994 Truss Set Screws	ME-6995 Road Bed Spares
ME-6996 Cord Lock Spares	ME-6997 Full Round (XYZ) Connector Spares
ME-6998 Axle Spares	ME-6999A Angle Connector Spares
PS-2198 Load Cell Amplifier	PS-2199 Load Cell and Amplifier Set
PS-2200 100 N Load Cell	PS-2201 5 N Load Cell
PS-2205 Dual Load Cell Amplifier	



