


Fan Rotation Base

ME-1251

Introduction

The Fan Rotation Base is designed to hold any fan compatible with PASCO carts and features a rotating platform, which can be used to demonstrate the vector nature of force. The base includes raised markings which indicate the angle the fan makes with the path of the cart. The accessory is designed to attach to any PASCO dynamics cart, such as the Smart Cart (ME-1240 or ME-1241) or the PAScar (ME-6933 or ME-6934)

Features

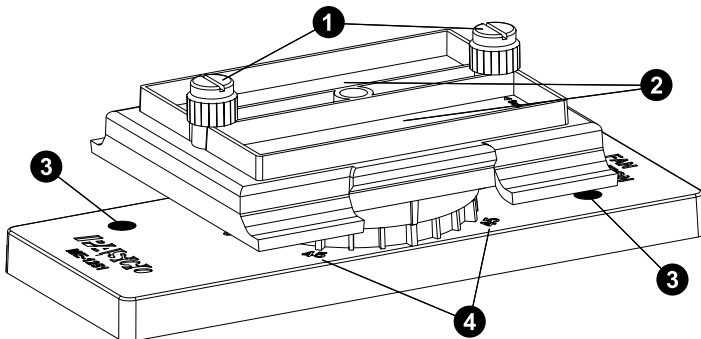

1 Mounting screw storage holes

Store the mounting screws here when they are not being used to secure the Fan Rotation Base to the cart.

2 Cart mass slots

Each of these slots can be used to hold a Compact Cart Mass (ME-6755), increasing the total mass of the cart and base system.

3 Mounting holes

Insert the mounting screws into these holes and turn clockwise to secure the base in place on a cart.

4 Numerical markings


Use to identify the angle of the fan with respect to the track. A raised line on the rotating disk indicates the current angle.

Required equipment:

- PASCO cart, such as the Smart Cart (ME-1240 or ME-1241) or the PAScar (ME-6933 or ME-6934)
- Any PASCO cart fan, such as the Smart Fan Accessory (ME-1242)

Usage

Mounting the apparatus

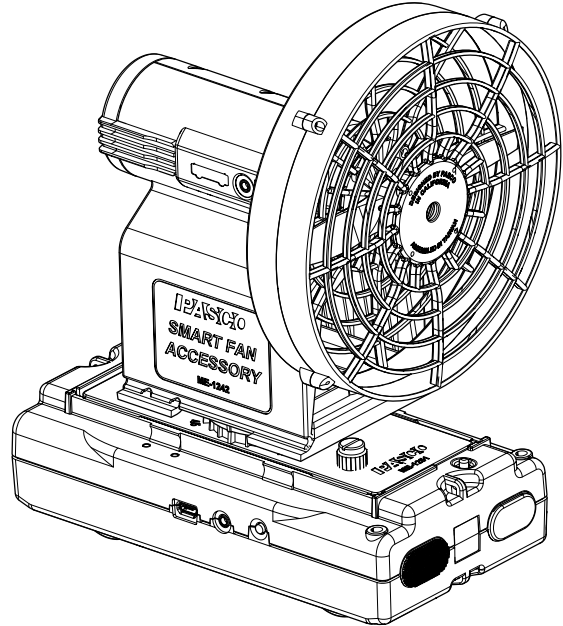

1. Remove the mounting screws from the storage holes on top of the rotating platform.
2. Place the Fan Rotation Base into the accessory tray of the cart.
3. Insert the mounting screws into the mounting holes. Tighten to secure the base in place.
4. Attach the fan to the attachment points on the rotating platform. A small amount of downward force may be required to secure the fan in place.

Rotating the base

The base is designed to rotate in 15 degree intervals, snapping between each position. The flat base features raised lines to indicate these positions, as well as numerical markings for 0° and 45°. A single raised line on the disk aligns with these markings, indicating the angle of the fan.

Technical support

Need more help? Our knowledgeable and friendly Technical Support staff is ready to answer your questions or walk you through any issues.

- Chat [pasco.com](https://www.pasco.com)
- Phone 1-800-772-8700 x1004 (USA)
+1 916 462 8384 (outside USA)
- Email support@pasco.com

Limited warranty

For a description of the product warranty, see the Warranty and Returns page at www.pasco.com/legal.

Copyright

This document is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Trademarks

PASCO and PASCO scientific are trademarks or registered trademarks of PASCO scientific, in the United States and in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.