




Goniometer

PS-2138, PS-2137

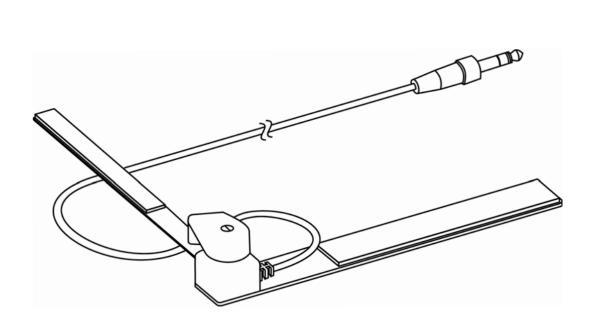

For Educational Use Only! Note: This is not a medical device. It is designed for educational use only and should not be used in any medical process such as life support or patient diagnosis. It is also not intended for use in graduate research or industry incluing industrial control or any type of industrial testing.

Goniometer PS-2138

Contents


Introduction
Probe Mounting
Sensor Setup
Suggested Activities
Experiment9
Experiment Teachers' Notes and Sample Data19
Other Sample Data
Safety
Specifications
Technical Support
Copyright
Limited Warranty



PS-2138 Goniometer

Goniometer

PS-2137, PS-2138

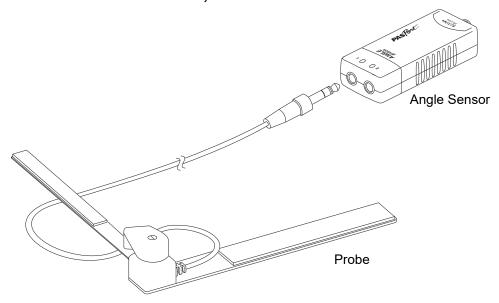

Equipment Included	Part Number	Part Number	
Probe	PS-2138		
Mounting straps (two small, one large)	PS-2547		
Angle Sensor (included in PS-2137 only)	PS-2139		

Model PS-2137 includes probe, straps and angle sensor; model PS-2138 includes probe and straps only.


Required or Optional Equipment

PASCO Interface and Data Collection Software	See PASCO catalog or www.pasco.com	
Second Goniometer Probe (optional)	PS-2138	

Introduction

With the PASPORT Goniometer students can analyze motions such as walking, running, throwing and kicking. They can also explore the physics of rotational dynamics using their arms and legs as experimental apparatus.

The Goniometer, in conjunction with a PASCO interface, measures and records the angle, angular velocity and angular acceleration of an elbow, knee or hip. The sensor can be used with a single Goniometer Probe (included) or with an optional second probe to measure two joints simultaneously. You can use multiple sensors together to support even more probes.

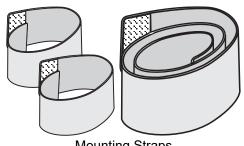


Mounting Straps

The probe is easily attached to the body using the included hook-and-loop mounting straps.

Goniometer Probe Mounting

The Goniometer Probe consists of two arms and a potentiometer. As the angle between the arms changes, the resistance of the potentiometer changes. The Angle Sensor, connected to the probe, measures the resistance of the potentiometer and converts it to an angle measurement. The sensor also calculates the angular velocity and angular acceleration from the rate at which the angle changes. The data is sent digitally to your PASCO interface at up to 100 samples per second.

The probe measures zero degrees (or radians) when it is fully open. A clockwise rotation of the narrow arm relative to the wide arm (as pictured) is measured as increasing angle.

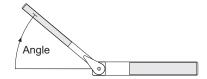

Probe Mounting

The mounting straps can be used in two ways. The easier method is to place the straps on the limbs, then stick the probe to the outside of the straps. For more secure attachment, tuck the probe arms inside the overlapping portions of the straps.

When mounting the Goniometer probe, the wide and narrow arms of the probe are interchangeable; the only difference will be the sign of the data collected. Attach the wide arm of the probe to the subject's upper arm when used on the left elbow, and to the subject's forearm when used on the right elbow. This will result in a flexion of either joint measured as a positive displacement. Similar reversals can be applied to knee and hip mounting.

Elbow

Place one strap around the upper arm midway between the elbow and shoulder. Place a second strap around the lower arm midway between the elbow and wrist.

With the hand relaxed, bend the elbow at a right angle. Align the probe's hinge with the elbow. Attach one arm of the probe to the subject's upper arm parallel to the humerus.



Right



Left

Extend the elbow and attach the other probe arm parallel to the ulna. Flex and extend the elbow a few times to check for proper alignment.

PS-2138 Sensor Setup

Knee

Place the large strap around the thigh just above the knee. Place a small strap around the upper part of the calf (combine both small straps end-to-end if necessary).

Bend the knee at a right angle. Align the probe's hinge with the knee. Attach one arm of the probe to the thigh parallel to the femur.

Extend the knee and attach the other probe arm parallel to the tibia.

Flex and extend the knee to check for proper alignment. Have the subject stand on both feet and make sure that the probe does not shift significantly.

Hip

Place the large strap around the waist. Combine both small straps end-to-end (if necessary) and place them around the upper thigh.

Have the subject stand and place his or her foot on a chair so that the thigh is horizontal. Align the probe's hinge with the hip joint. Attach one arm of the probe to the thigh parallel to the tibia.

Have the subject stand on both feet; attach the other probe arm vertically to the waist strap.

Move the hip joint through its full range of forward and backward rotation to check for proper alignment.



Sensor Setup

Connect one or two Goniometer Probes to the Angle Sensor. Connect the Angle Sensor to a PASCO interface. The interface will collect data for angle, angular velocity, and angular acceleration from each probe. You can select units of degrees or radians in the software.

The default sampling rate of the sensor is 20 Hz. In most cases this is sufficient, but for faster movements you may wish to increase the rate to 50 or 100 Hz.

Optional Data Smoothing

The sensor calculates the angular velocity and angular acceleration from the measured angle data. The angular velocity is the change in angle between consecutive samples divided by the time between samples. Angular acceleration is the change between consecutive velocity calculations divided by the time between samples.

The Goniometer is very sensitive to small variations in the angular velocity, so you may see a lot of variation in angular acceleration. To make the angular acceleration data easier for students to interpret, use the smooth function in the PASCO data collection software. See the Online Help for the software for more information.

Optional Calibration

The Goniometer does not normally require calibration. To increase the accuracy of measurements made over a limited range of motion, the Goniometer may be manually calibrated. In the data collection software, click the Calibrate button in the t Setup window. Set the probe at a known angle, enter the angle under Point 1 and click the Set button. Set the probe at another known angle, enter the angle under Point 2 and click the Set button.

Suggested Activities

Analysis of Gait and other motions

Collect angle data of the knee while walking. Does it approximate simple harmonic motion? Explain what you observe.

Collect angle data of the hips during walking, fast walking and running.

- How does the angle of forward rotation compare to the angle of backward rotation?
- Compare the left and right hips. Are they symmetrical?
- Compare the range of movement and period of oscillation for walking, fast walking and running? What patterns do you observe?
- Compare data from different students walking at the same speed (walk sideby-side or use a motion sensor to monitor speed).
- For that class, make histograms of range of motion and period of oscillation. Do any patterns emerge?
- Make graphs of range of motion and period vs. height. Is there a correlation?

Measure the period of the leg swinging freely and compare it to the period of oscillation when the subject is walking at his or her most comfortable pace. For the class, investigate the relationship between leg length, L, and walking period, \mathcal{T} . For all pendulums $T \propto \sqrt{L}$; is this true for human legs*?

Analyze non-periodic movements such as throwing, kicking, and lifting. When performing the movements, move only the joint that is being measured. You can collect data on the linear motion of lifted, kicked and thrown objects using Photogate Tape (ME-6664), a Photogate (ME-9204B) and a Digital Adapter (PS-2159).

Collect data from two or more joints simultaneously while walking, running, jumping, throwing, kicking, etc. How do the joints work together when performing these actions?

*For a detailed analysis of the leg as a physical pendulum see: A. Dumont and C. Waltham, 1997, Walking, *The Physics Teacher*, 35 (6): 372– 376

Oscillations

Collect data for the following:

- Lower leg freely dangling about the knee showing simple harmonic motion, subject seated on a high surface.
- Leg with unbent knee freely dangling about the hip, subject standing on the opposite foot on a low surface.
- · Leg with knee bent at right angle freely dangling about the hip.

Do angle, angular velocity, and angular acceleration approximate simple harmonic motion? Determine the period, frequency and amplitude of the oscillations.

What is the relationship between the phases of angle, angular velocity, and angular acceleration?

How does bending the knee affect the frequency of the dangling leg?

Skeletal Parts

Mount the Goniometer on articulated skeletal parts. In conjunction with force sensors attached with string the points of tendon attachment, measure the forces exerted by muscles when lifting objects of various mass, or performing throwing and kicking movements.

PS-2138 Experiment

Experiment

In this activity you will rotate your arm about your elbow and investigate how the position, velocity and acceleration of your hand relate to the angle, angular velocity and angular acceleration of your arm.


Equipment	Part Number
Goniometer Probe	PS-2138
Angle Sensor	PS-2139
PASCO interface and Data Collection Software	See PASCO catalog or www.pasco.com
Motion Sensor	PS-2103
Acceleration Sensor	PS-2118, PS-2119 or PS-2136
Wall-mounted white board and pen	

Part 1: Arc Length vs. Angle

When you rotate your forearm about your elbow (while keeping your upper arm stationary) your hand traces out an arc, or part of a circle. You will use a pen and white board to mark the path of your hand. The Goniometer will measure the angle through which your forearm rotates.

Setup

- · Connect the Goniometer to the interface.
- Mount the Goniometer on your elbow* so that a flexion of the joint is measured as a positive angular displacement (as pictured below).

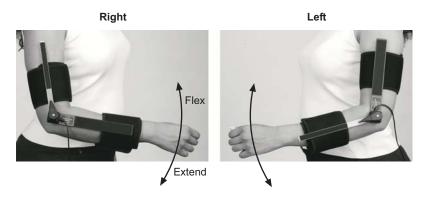

*If you are right-handed, mount the Goniometer on your right elbow; if you are left-handed, mount it on your left elbow.


1. Stand next to the white board with your arm relaxed at your side. Let your elbow and the *back* of your hand touch the board.

Goniometer Experiment


2. Place a pen in your hand (as shown) so that you will draw on the board while keeping the back of your hand closest to the board.

Have a partner mark the location of your elbow on the board, and measure the distance from your elbow to the pen.

4. Distance from elbow to pen

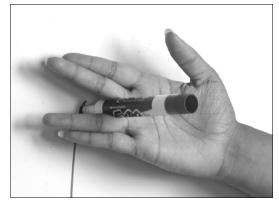

It is important not to twist your hand in order to allow your arm to rotate in a plane parallel to the board

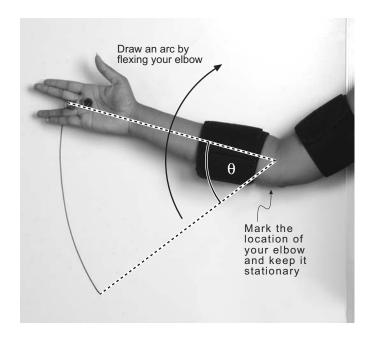

Procedure

- 1. Fully extend your elbow and place the pen tip on the board.
- 2. Start data collection.
- **3.** Draw an arc on the board by flexing your elbow. Move only your forearm and hand, while keeping your elbow at the marked location on the board.
- **4.** Stop data collection.

Analysis

1.	Look at the graph of Angle vs. Time . According to the graph, what angle ir
	radians did you trace out on the board?

2. When measured in radians, the arc angle (θ) is the ratio of the arc length (s) to radius (r). In this case r is the distance from your elbow to the pen.

$$\theta = s/r$$


According to this theoretical relationship, and your measured values of θ and r, how far did your hand travel?

$$s =$$
_____(theoretical)

3. Measure the length of the arc that you drew and record it below. (Tape a piece of string to the board exactly over the arc. Mark the endpoints of the arc on the string, then un-tape the string, lay it straight on a table, and measure the distance between the marks.)

$$s =$$
_____ (actual)

4.	How does the theoretical value of s compare to the actual distance that your
	hand traveled?

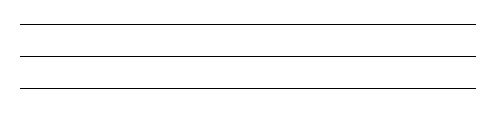

Goniometer Experiment

Part 2: Tangential Velocity vs. Angular Velocity

When you rotate your forearm about the elbow, your hand does not move in a straight line, but it always moves in a direction perpendicular to your forearm. This direction is described as tangential.

As you rotate your arm, the magnitude of your hand's tangential velocity (v_{τ}) equals the change in the arc length traced by your hand (Δs) divided by the change in time (Δt)

$$v_{\tau} = \Delta s / \Delta t$$

The angular velocity (ω) of your forearm equals the change in angle ($\Delta\theta$) divided by the change in time (Δt).

$$\omega = \Delta\theta/\Delta t$$

In Part 1, you discovered that the relationship between arc length and angle is $s = r \theta$, thus:

$$\Delta s = r \Delta \theta$$

Predict

the ab	ou will use the Mo e Goniometer to r ove equations, w ationship betwee	neasure the ang rite an equation	ular velocity of	your forearm. B	ased on the	and
_						
_						


Setup

- 1. Connect the Goniometer and Motion Sensor to the interface (or interfaces).
- 2. Place the Goniometer on your elbow so that a flexion of the joint is measured as a positive angular displacement.

PS-2138 Experiment

3. Mount the Motion Sensor in front of you at shoulder level and about 60 cm from your chest.

4. Hold your arm (as shown in the picture) so that your forearm will rotate in a horizontal plane with your hand in front of the Motion Sensor.

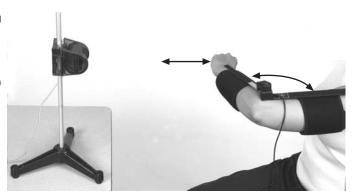

Procedure

- 1. Start data collection.
- 2. While keeping your shoulder still, bend your elbow to move your hand toward the Motion Sensor, then move your hand back toward your chest. (Don't let your hand get closer than 15 cm from the sensor.)
- 3. Stop data collection.

Analysis

In this analysis, use units of m/s for velocity and rad/s for angular velocity.

1.	Look at graphs of Velocity vs. Time and Angular Velocity vs. Time together. How are the two graphs related?	_
		- - -
2.	Create a graph of Velocity vs. Angular Velocity .* Qualitatively describe the graph.	*Plot Velocity on the vertical axis and Angular Velocity on the horizontal axis.
3.	Apply a linear fit to the graph. What is the slope of the best-fit line? Include units.* $Slope = ___$	*When determining the units, remember to consider the definition of a radian. Though a radian is usually used as a unit of measure, it is actually a unitless quantity. Thus m/rad is equivalent to m.

4.	Write the equation of the best-fit line in terms of $v_{_T},\omega$ and slope. Does the relationship represented by this equation support the prediction that you made earlier?
5.	What physical quantity is represented by the slope? Measure this quantity directly and compare it to the value of slope.

PS-2138 Experiment

Part 3: Centripetal and Tangential Accelerations

When your forearm rotates about your elbow the velocity of your hand is entirely in the tangential direction, but the acceleration is not. There are two components to the acceleration of your hand: tangential (perpendicular to your forearm), and centripetal (parallel to your forearm.) You will use the Acceleration Sensor to measure both components.

The magnitude of tangential acceleration (a_{τ}) is equal to the change in magnitude of tangential velocity (Δv_{τ}) divided by the change in time (Δt) .

$$a_{\scriptscriptstyle T} = \Delta v_{\scriptscriptstyle T} / \Delta t$$

The angular acceleration (α) of your forearm equals the change in angular velocity ($\Delta\omega$) divided by the change in time (Δt).

$$\alpha = \Delta \omega / \Delta t$$

In Part 2, you discovered that the relationship between tangential velocity and angular velocity is $v_{\tau} = r \omega$, thus:

$$\Delta v_{T} = r \Delta \omega$$

Predict


You will use the Acceleration Sensor to measure the tangential acceleration of your hand, and the Goniometer to measure the angular acceleration of your forearm. Based on the above equations, write an equation predicting what you will discover about the relationship between $a_{\scriptscriptstyle T}$ and α .

Setup

- 1. Connect the Goniometer and Acceleration Sensor to the interface (or interfaces).
- 2. Set the sample rates of both sensors to 20 Hz.
- **3.** Place the Goniometer on your elbow so that a flexion of the joint is measured as a positive angular displacement.

Goniometer Experiment



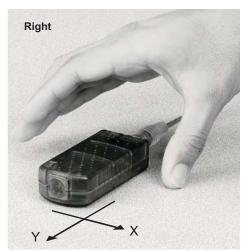



- 4. Hold the Acceleration Sensor in your hand with the sensor's X-axis parallel to your forearm and pointing toward your elbow, and the Y-axis orthogonal to your forearm and pointing in the direction that your hand moves when you flex your elbow (as shown in the picture).
- **5.** Measure the distance, *r*, from your elbow to the Acceleration Sensor.

r =

6. Hold your arm so that your forearm will rotate in a horizontal plane (as pictured), and the X-Y plane of the Acceleration Sensor is also horizontal.

Procedure

- **1.** Extend your elbow.
- 2. Start data collection.
- 3. Quickly flex your elbow, wait a moment, then quickly extend your elbow.*
- 4. Stop data collection.

*The Acceleration Sensor must remain level (its X-Y plane parallel to the floor) throughout data collection. PS-2138 Experiment

Analysis

In this analysis, use units of m/s/s or m/s^2 for acceleration, rad/s for angular velocity, and rad/s/s or rad/s^2 for angular acceleration.

Tangential Acceleration

1.	Look at graphs of Angular Acceleration vs. Time and Tangential Acceleration vs. Time together.* How are the graphs related?	*Tangential Acceleration is measured by the sensor as "Acceleration, Y."
		-
2.	Create a graph of Tangential Acceleration vs. Angular Acceleration and apply a linear fit.* What is the slope of the best-fit line (including units)?	*Plot Tangential Acceleration on the vertical axis and Angula
3.	Write the equation of the best fit line in terms of $a_{\rm T}$, α and slope. Does the relationship represented by the best-fit line support the prediction that you made earlier?	Acceleration on the horizontal axis.
		-
4.	What physical quantity is represented by the slope? Measure this quantity directly and compare it to the value of slope.	_
		- - -

Goniometer Experiment


Centripetal Acceleration

•	Look at graphs of Angular Velocity vs. Time and Centripetal Acceleration vs. Time together.* How are the graphs related?	*Centripetal Acceleration is measured by the sensor as "Acceleration, .X."
)	The Acceleration Sensor was oriented so that acceleration toward the elbow was measured as positive. When the angular velocity of your arm was positive, was the centripetal acceleration of your hand toward or away from	-
3.	when the angular velocity of your arm was <i>negative</i> , was the centripetal acceleration of your hand toward or away from your elbow?	
1.	Create a graph of Centripetal Acceleration vs. Angular Velocity .* The theoretical relationship between Centripetal Acceleration (a_c) and Angular Velocity (ω) is $a_c = r \omega^2$, where r is a constant. Does your data appear to support this relationship? Explain your reasoning.	*Plot Centripetal Acceleration on the vertical axis and Angular Velocity on the horizonta axis.
5.	Create a graph of a_c versus ω^2 . Apply a linear fit. According to the slope of the best-fit line, what is the value of r (including units)?	_
3 .	$r = \underline{\hspace{1cm}} \text{ (from best-fit line)}$ Compare this value to the distance from your elbow to the acceleration sensor.	
		-
		_

18

Experiment Teachers' Notes and Sample Data

Part 1

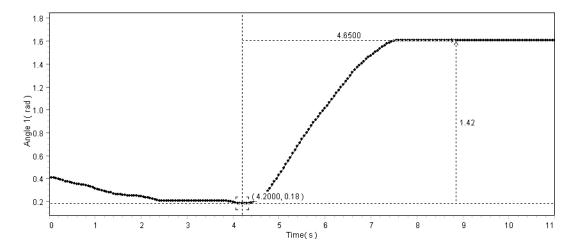


In this example r = 0.38 m.

- **1.** Arc Angle, θ = 1.42 rad
- **2.** $s = r \theta = (0.38 \text{ m}) (1.42 \text{ rad}) = 0.54 \text{ m}$, theoretical
- **3.** s = 0.57 m, actual
- **4.** In this example the theoretical and actual values of s differ by 5%.


Part 2

Students should predict $v_{\tau} = r_{\odot}$

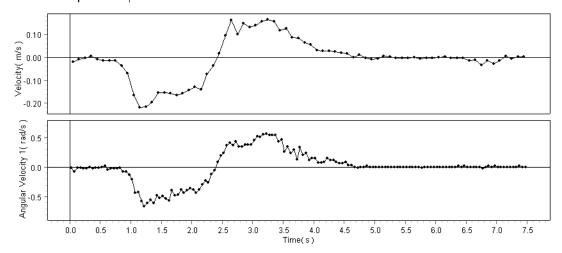

19

1. The graphs of Velocity vs. Time and Angular Velocity vs. Time appear to be directly proportional.

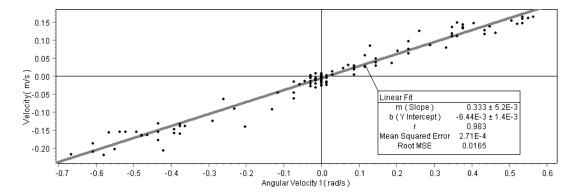


- **2.** The graph of Velocity vs. Angular Velocity shows a directly proportional, linear relationship.
- 3. In this example slope = $0.333 \text{ m} \pm 0.005 \text{ m}$
- **4.** v_{τ} = slope × ω v_{τ} = (0.333 m) ω This equation supports the prediction.
- **5.** Slope is *r*, the distance from the elbow to the hand. In this case the actual value of *r* is 0.36 m. The theoretical value (from slope) and actual value (from direct measurement) differ by about 8%.


Note that the directly measured values of r in Part 1 and Part 2 are slightly different, though they are from the same student. This is due to the different hand positions used in each part.

Part 3


Students should predict $a_{\tau} = r\alpha$

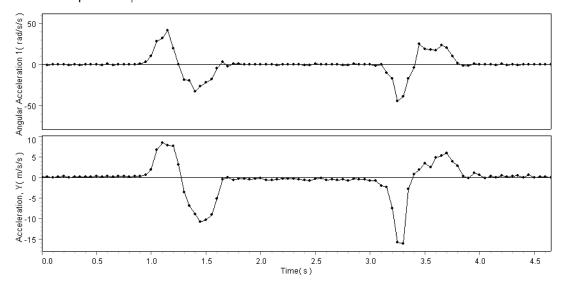


1. The graphs of Angular Acceleration vs. Time and Tangential Acceleration vs. Time appear to be directly proportional.

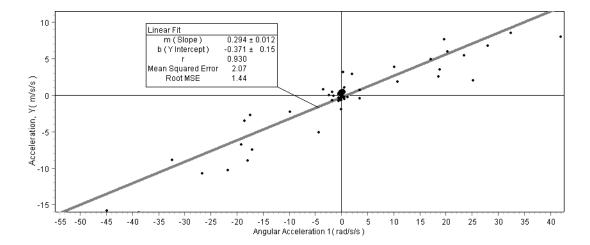


- 2. In this example, the slope of the best fit line is $0.29 \text{ m} \pm 0.01 \text{ m}$.
- 3. $a_{\scriptscriptstyle T}$ = slope × α $a_{\scriptscriptstyle T}$ = (0.29 m) α This equation supports the prediction.
- **4.** Slope is equal to *r*. In this case the actual value is 0.36 m. The theoretical value (from slope) and actual value (from direct measurement) differ by about 24%.

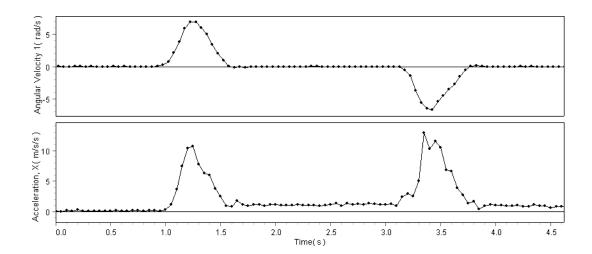

- **5.** The graphs of Angular Velocity vs. Time and Centripetal Acceleration vs. Time appear to show a proportionality between a_c and the *magnitude* of ω .
- **6.** When ω was positive, a_c was also positive, therefor the centripetal acceleration was toward the elbow.
- 7. When ω was negative, $a_{\rm c}$ was positive, therefor the centripetal acceleration was toward the elbow.

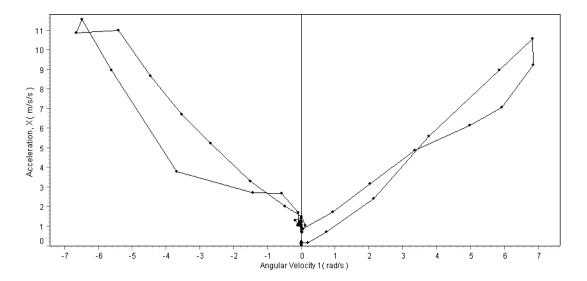

8. In theory, the graph of a_c versus ω is a parabola. Though the data contains a lot of scatter, that relationship is evident in the collected data.

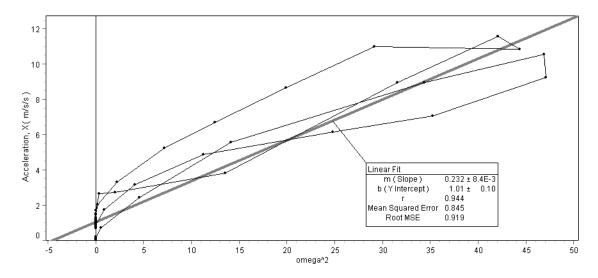


- 9. On the graph of a_c versus ω^2 , the slope of the best-fit line is the theoretical value of r, in this case 0.232 m \pm 0.008 m.
- **10.** The theoretical and actual values of *r* differ by about 55%.


In Part 3, your students will probably find that the values of *r* according to the best-fit lines differ significantly from the actual value; and that the graphs themselves contain significant scatter or noise.

Have them compare the different sensors, procedures, and mathematics (including the calculations done inside the sensor) used in each part of the experiment, and consider how these factors contribute to precision and accuracy of the collected data.


Other Sample Data

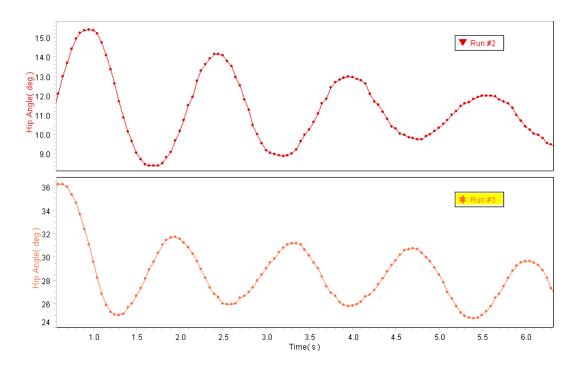


Simple Harmonic Motion of the leg rotating about the hip, dangling freely and oscillating at its natural frequency, with the knee unbent (top) and bent (bottom). Note the higher frequency of the leg with bent knee.

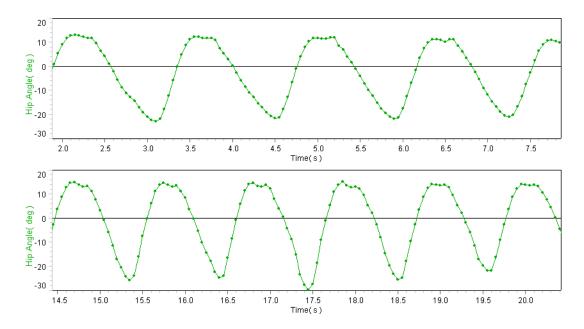

Hip angle while waking at a normal pace (top) and walking quickly (bottom). Note the difference in frequency *and* amplitude.

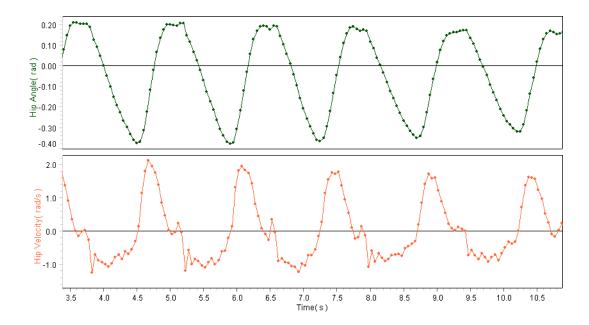

The angle (top) and angular velocity (bottom) of the hip while walking. The maximum slope of the angle plot is about 1.9 rad/s, equal to the maximum value of angular velocity.

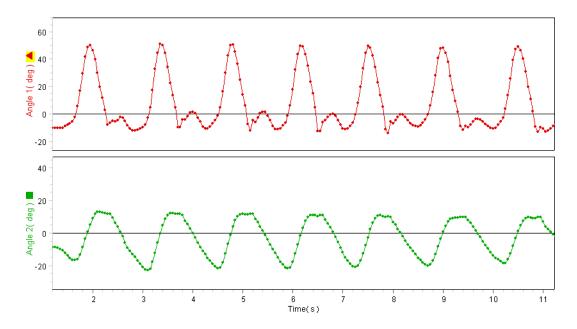

Angle data from the knee (top) and hip (bottom) of a walking subject, acquired using two probes simultaneously. Note the phase relationship between the joints.

PS-2138 Goniometer

Safety

Read the instructions before using this product. Students should be supervised by their instructors. When using this product, follow the instructions in this manual and all local safety guidelines that apply to you.

Ensure that students are aware of these safety precautions before using the Goniometer.

Place the mounting straps on the body snugly but not too tightly. The straps should not constrict blood flow or breathing. Remove all straps at the first sign of discomfort.

When performing physical activities such as walking or running, do so only at a comfortable rate. Some activities call for students to stand or sit on raised surfaces; use only surfaces that are safe and appropriate for these activities.

Specifications

PS-2138 Goniometer Probe used with PS-2139 Angle Sensor and a PASPORT interface.

Range	±170°			
Accuracy	±1° calibrated ±3° uncalibrated			
Resolution	0.042°			
Sampling rate	20 Hz default 500 Hz maximum			
Probe arm length	21 cm			
Mounting straps	Large: 15 × 120 cm Small: 10 × 18 cm			

Author: Alec Ogston

The PASCO Goniometer was developed in cooperation with Dr. Nancy Beverly, Mercy College, Dobbs Ferry NY.

Technical Support

For assistance with any PASCO product, contact PASCO at:

Address: PASCO scientific

10101 Foothills Blvd.

Roseville, CA 95747-7100

Phone: (916) 462-8384

(800) 772-8700

Web: www.pasco.com
Email: techsupp@pasco.com

Limited WarrantyFor a description of the product warranty, see the PASCO catalog.

Copyright

The PASCO scientific 012-08904C *Goniometer Instruction Manual* is copyrighted with all rights reserved. Permission is granted to non-profit educational institutions for reproduction of any part of this manual, providing the reproductions are used only in their laboratories and classrooms, and are not sold for profit. Reproduction under any other circumstances, without the written consent of PASCO scientific, is prohibited.

Trademarks

PASCO, PASCO scientific, and PASPORT are trademarks or registered trademarks of PASCO scientific, in the United States and/or in other countries. All other brands, products, or service names are or may be trademarks or service marks of, and are used to identify, products or services of, their respective owners. For more information visit www.pasco.com/legal.

Product End of Life Disposal Instructions:

This electronic product is subject to disposal and recycling regulations that vary by country and region. It is your responsibility to recycle your electronic equipment per your local environmental laws and regulations to ensure that it will be recycled in a manner that protects human health and the environment. To find out where you can drop off your waste equipment for recycling, please contact your local waste recycle/disposal service, or the place where you purchased the product.

The European Union WEEE (Waste Electronic and Electrical Equipment) symbol (to the right) and on the product or its packaging indicates that this product must not be disposed of in a standard waste container.

