Instruction Sheet for the PASCO Model WA-9870

Wave Driver Post (2) Installation Instructions

Introduction

The WA-9870 Wave Driver Post provides a set of two posts to be glued to any 5 1/4-inch speakers as indicated in Figure 1.

Equipment required but not included:

- Glue; preferrably RTV (black or clear)

Assembling the Plumb Bob

- 1. Carefully draw an even bead of RTV around the internal perimeter of the speaker.
- 2. Allow the RTV to dry according to manufacturer's recommendations before setting the post in place. Assure that the seal around the post is complete.

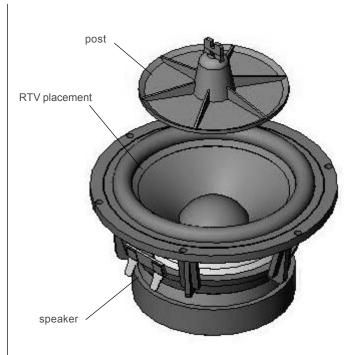


Figure 1. Alignment of Post on Speaker

© 2004 PASCO scientific



10101 Foothills Blvd. • P.O. Box 619011 • Roseville, CA 95678-9011 USA Phone (916) 786-3800 • FAX (916) 786-8905 • web: www.pasco.com

